IKA

designed for scientists

Betriebsanleitung Ursprungssprache	DE	3
Operating instructions	EN	6
Mode d'emploi	FR	9
Руководство по эксплуатации	RU	12

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Contents

	Page
Explanation of symbols	6
Safety instructions	6
Correct use	7
Unpacking	7
Useful information	7
Commissioning	7
Maintenance and cleaning	8
Warranty	8
Technical data	8

Explanation of symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the instrument or to system components.

This symbol indicates information **which is important for ensuring that the operations of the instrument are performed efficiently and for using the instrument**. Failure to observe this information can result in inaccurate results.

Danger - Risk of hand injury.

Safety instructions

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Follow the safety instructions, guidelines, occupational health, safety and accident prevention regulations.
- Set up the lab lift in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- Protect the instrument from bumps and impacts.
- The feet of the appliance must be clean and undamaged.
- Position one laboratory device or vessel in the centre and several devices or vessels evenly on the working plate in use.

Never put overweight load on the lift, see section "**Technical data**" for the maximum load.

- Pay attention to the device or vessel on the working plate when increase or decrease the height of the lab lift.

Pay attention to hand injury when operate the lab lift.

Correct use

• Use

For raising / supporting laboratory devices or vessels.

Intended use: Tabletop device.

• Range of use

- Laboratories
- Pharmacies
- Schools
- Universities

This instrument is suitable for use in all areas except industry areas.

The safety of the user cannot be guaranteed if the instrument is operated improperly or contrary to the manufacturer's specifications.

Unpacking

• Unpacking

- Please unpack the device carefully
- In the case of any damage a fact report must be sent immediately (post, rail or forwarder).

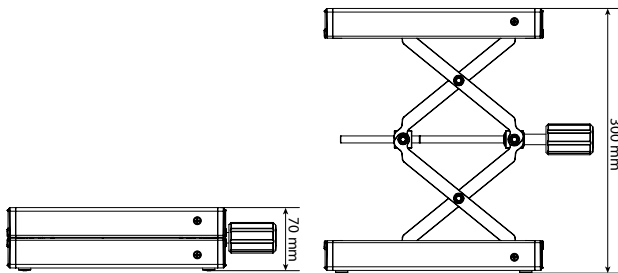
• Delivery scope

- IKA® LABLIFT m
- An operation instructions
- A warranty card
- A decontamination clearance certificate.

Useful information

IKA® LABLIFT m is suitable for raising and supporting small laboratory devices or vessels. The maximum raising weight can reach up to 7 kg (**Max. dynamic load**) and supporting weight up to 25 kg (**Max. static load**). The working height can reach from 70 mm to 300 mm.

The handy shape handwheel ensure the operation easily and comfortably. The self-lubricating bearing reduces the friction largely. The auxiliary springs ensure the user could raise large load easily and smoothly especially in the lowest position.

Commissioning

Place the IKA® LABLIFT m on working table firmly. Turn the handwheel in clockwise and counterclockwise direction, the working plate of the device goes up and down accordingly.

Observe the "**Max. dynamic load**" and "**Max. static load**" listed under "**Technical data**".

Maintenance and cleaning

The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

Only use cleansing agents which have been recommended by **IKA®**.

Dirt

Dye

Construction material

Cosmetics

Foodstuffs

Fuel

Cleaning agent

isopropyl alcohol

water containing tenside /

isopropyl alcohol

water containing tenside /

isopropyl alcohol

water containing tenside

water containing tenside

For materials which are not listed, please request information from **IKA®** application support.

Wear protective gloves during cleaning the instruments.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

Spare parts order

When ordering spare parts, please give:

- machine type
- manufacturing number, see type plate
- item and designation of the spare part, see **www.ika.com**, spare parts diagram and spare parts list.

Repair

Please send in instrument for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For this you should request the "Decontamination Clearance Certificate" from **IKA®**, or use the download printout of it from the **IKA®** website **www.ika.com**.

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited