
# IKA

designed for scientists

## IKA HG 600 IKA HG 1000

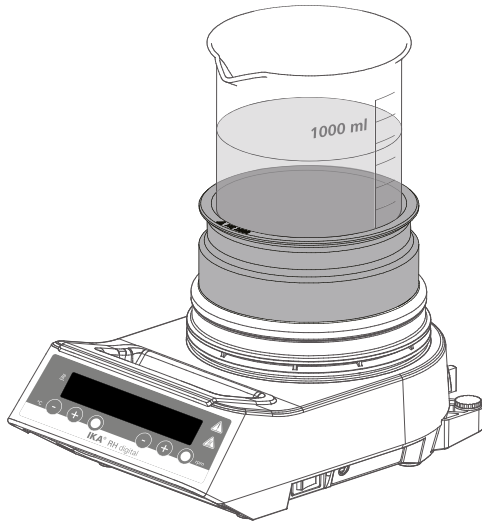Technische Information  
Technical information

DE  
EN

Ursprungssprache

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## Explication of warning symbols

-  **DANGER** Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.
-  **WARNING** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
-  **CAUTION** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
-  **NOTICE** Indicates practices which, if not avoided, can result in equipment damage.
-  **WARNING** Reference to exposure to a hot surface!
-  Hot surface!

## Safety instructions

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

 **WARNING** During handling the heating jackets, hot surface can cause risk of burning! Wear gloves when handling the heating jackets!

- Pay attention to the residual heat after processing.
- Never move the heating jackets as long as they are hot.
- Observe the operating instructions of the respective magnetic stirrer.

 **NOTICE** Beware of hazards due to:

- flammable materials
- combustible media with a low boiling temperature
- glass breakage
- unsafe condition of container.

- The heating jacket and the beaker must be properly placed on the heating plate and must not come off by themselves. The center of gravity of the heating jacket and the beaker must lie in the heating plate.

## Proper use

**IKA HG 600** or **HG 1000** heating jacket can be used for matching with **IKA** magnetic stirrers with heating or heating plate. The heating jacket improves heat transfer into the medium. It can cut down the heating time considerably.

Anodizing color of the heating jacket will change with aging, it do not affect normally heating function of the heating jacket.

## Cleaning

 **WARNING** Pay attention to the residual heat prior to cleaning. Never clean the devices when they are still hot!

Use only cleaning agents which have been approved by **IKA** to clean the device: water containing tenside / isopropyl alcohol. Wear protective gloves while cleaning the device. If a different cleaning or decontamination method than the method defined by **IKA** is planned, the user must ascertain with **IKA** that this method does not damage the device.

**Note:** Please don't clean the devices in a lab dishwasher! Or else, the coating may peel off, but has not affect to the functionality of the devices.

## Technical data

		HG 600	HG 1000
For beaker volume	ml	600	1000
Material		Aluminium	
Inner diameter	mm	91	106
Dimensions	mm	Ø 111 x 50	Ø 126 x 50
Weight	kg	0.35	0.40

*Subject to technical changes!*

## Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs. The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

# IKA

designed for scientists

**IKA-Werke GmbH & Co. KG**  
Janke & Kunkel-Straße 10, 79219 Staufen, Germany  
Phone: +49 7633 831-0, Fax: +49 7633 831-98  
eMail: sales@ika.de

**USA**  
**IKA Works, Inc.**  
Phone: +1 910 452-7059  
eMail: usa@ika.net

**KOREA**  
**IKA Korea Ltd.**  
Phone: +82 2 2136 6800  
eMail: sales-lab@ika.kr

**BRAZIL**  
**IKA Brazil**  
Phone: +55 19 3772 9600  
eMail: sales@ika.net.br

**MALAYSIA**  
**IKA Works (Asia) Sdn Bhd**  
Phone: +60 3 6099-5666  
eMail: sales.lab@ika.my

**CHINA**  
**IKA Works Guangzhou**  
Phone: +86 20 8222 6771  
eMail: info@ika.cn

**POLAND**  
**IKA Poland Sp. z o.o.**  
Phone: +48 22 201 99 79  
eMail: sales.poland@ika.com

**JAPAN**  
**IKA Japan K.K.**  
Phone: +81 6 6730 6781  
eMail: info\_japan@ika.ne.jp

**INDIA**  
**IKA India Private Limited**  
Phone: +91 80 26253 900  
eMail: info@ika.in

**UNITED KINGDOM**  
**IKA England LTD.**  
Phone: +44 1865 986 162  
eMail: sales.England@ika.com

**VIETNAM**  
**IKA Vietnam Company Limited**  
Phone: +84 28 38202142  
eMail: sales.lab-vietnam@ika.com

Discover and order the fascinating products of IKA online:  
[www.ika.com](http://www.ika.com)

 IKAworldwide  IKAworldwide /// #lookattheblue  @IKAworldwide