IKA

designed for scientists

IKA lab disc

Betriebsanleitung	DE	4
Operating instructions	EN	8
Mode d'emploi	FR	12
使用说明	ZH	16
Indicaciones de seguridad	ES	20
Veiligheidsinstructies	NL	21
Normi di sicurezza	IT	23
Säkerhetsanvisningar	SV	24
Sikkerhedshenvisninger	DA	26
Sikkerhedsanvisninger	NO	27
Turvallisuusohjeet	FI	29
Normas de segurança	PT	30
Wskazówki bezpieczeństwa	PL	32
Bepečnostní upozornění	CS	33
Biztonsági utasítások	HU	35
Varnostna navodila	SL	36
Bezpečnostné pokyny	SK	38
Ohutusjuhised	ET	39
Drošības norādes	LV	41
Saugos reikalavima	LT	42

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Contents

	Page
Safety instructions	8
Operator's control	9
Unpacking	9
Intended use	10
Starting up	10
Maintenance and cleaning	10
Accessories	11
Warranty	11
Technical data	11

Safety instructions

For your protection

- **Read the operating instructions in full before starting up and follow the safety instructions.**

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Caution - Magnetism! Effects of the magnetic field have to be taken into account (e.g. data storage media, cardiac pacemakers...).

- Wear your personal protective equipment in accordance with the hazard category of the media to be processed. Otherwise there is a risk from:
 - splashing and evaporation of liquids
 - ejection of parts
 - release of toxic or combustible gases.

- Use only the AC adapter supplied.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The base plate must be clean and intact.
- Check the appliance and accessories beforehand for damage. Do not use damaged components.
- Reduce the speed if
 - the medium splashes out of the vessel because the speed is too high
 - the appliance is not running smoothly
 - the container moves on the base plate.
- Beware of and avoid hazards due to
 - flammable materials
 - combustible media with a low boiling temperature
 - incorrect container size
 - overfilling of media
 - unsafe condition of container.
- The appliance may heat up when in use.
- Process pathogenic materials only in closed vessels under a suitable extractor hood.
- Do not use the device in explosive atmospheres, it is not EX-protected.
- With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.
- To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any energy produced in other ways, e.g. through light irradiation.
- Safe operation is guaranteed only with the use of original IKA

accessories.

- The appliance must only be disconnected from the mains supply by pulling out the plug-in power supply unit.
- The socket for the the plug-in power supply unit must be easily accessible.
- The device will automatically restart following any interruption to the power supply.
- It may be possible for wear debris from rotating accessory parts to reach the material being processed.
- When using the PTFE-coated magnetic bars, the following has to be noted: Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline-earth metals, as well as with fineparticeld powders of metals of the 2. and 3. group of the periodical system at temperatures above 300-400 °C. Only elementary fluorine, chlorine trifluoride and alkaline metals do attack PTFE, halogen hydrocarbons have a reversible swelling effect. (Source: Römpps Chemie-Lexikon and „Ullmann“ Bd.19)

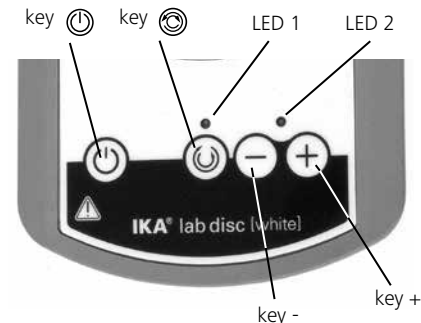
To the protection of the equipment

Caution: Equipment may not be poured over with aggressive media! Equipment never under water set!

- The voltage stated on the type plate must correspond to the mains voltage.
- Do not cover the device, even partially e.g. with metallic plates or film. This results in overheating.
- Protect the appliance and accessories from bumps and impacts.
- Ensure that the base plate is kept clean.

Operator's control

Unpacking

Unpacking

- Please unpack the device carefully
- In the case of any damage a detailed report must be sent immediately (post, rail or forwarder).

Scope of delivery

- Magnetic stirrer
- Power supply unit with exchangeable adapters
- Operating instructions

Intended use

• Use

- for mixing liquids

• Area of use

Indoor environments similar to that a laboratory of research, teaching, trade or industry area.

Starting up

- Connect the appliance to a socket (the voltage stated on the nameplate must correspond to the mains voltage).
- Number of revolutions to minimum value set (switch on with , with  reduce number of revolutions to LED 2 permanently shines)
- Place a magnetic rod in a vessel which is penetrable for magnetic field lines and fill the vessel with the medium to be processed.
- Place the vessel on the centre of the set-up surface.
- Set the speed using  and  (LED 2 will flash)

To elect the reverse mode and invert the direction of stirring at regular intervals, press the button . In the revers mode the stirring direction will be inverted every 30 seconds; there will be pause of 5 seconds between each changeover. The LED 1 will flash when the revers mode is active.

The speed setting and revers mode selection will remain unchanged when the instrument is unplugged.

Caution: At some speeds, irregularities in the shape of the stirring bar and of the bottom of the container can cause the bar to move irregularly during stirring. Increase or decrease the speed until the stirring bar begins to rotate evenly.

Maintenance and cleaning

The device is maintenance-free.

Cleaning

- For cleaning disconnect the mains plug!
- Use only cleaning agents which have been approved by IKA to clean IKA devices. These are water (with tenside) and isopropanol.
- Wear protective gloves during cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Accessories

IKA - stirring bars:

- TRIKA® 25
- IKAFLON® 30
- Stirring bar remover

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine directly to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

Technical data

Plug-in power supply unit:

Rated voltage	VAC	100 - 240
Rated frequency	Hz	50/60
Output voltage	VDC	12
Absorbed power	W	5
Output power	W	min. 3

Magnetic stirrer:

Speed range	rpm	15 - 1500
Stirred quantity (max.)	ml	800 (at 1500 rpm with 25mm stirring bar)

Operating time	%	100
Protection class		II
Pollution degree		2
Overvoltage category		II
EN 60529 protection class		IP 65
Ambient humidity (relative)	%	80
Ambient temperature	°C	5 to 40 (in the operation)
Transport temperature	°C	-25 to +70
Storage temperature	°C	-10 to +70
Material: Housing		PA GF

	TPU	
Front membrane sheet		Polyester
Weight	kg	0.3
Installed weight (max.)	kg	2.0
Set-up surface	mm	ø 100
Dimensions (W x D x H)	mm	116 x 175 x 12

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IKA Vietnam Company Limited
IKA India Private Limited