2000007192

WiCo LR 1000 control\_RET control-visc\_122015

# IKA®

## IKA® WiCo LR 1000 control

## IKA® WiCo RET control-visc

Betriebsanleitung	DE	02
Ursprungssprache		
Operating instructions	EN	10
Mode d'emploi	FR	18
Руководство пользователя	RU	26
使用说明	ZH	34
取扱説明書	JA	41
사용 설명서	KO	48
Indicaciones de seguridad	ES	55
Veiligheidsinstructies	NL	56
Norme di sicurezza	IT	58
Säkerhetsanvisningar	SV	59
Sikkerhedshenvisninger	DA	61
Sikkerhetsanvisninger	NO	62
Turvallisuusohjeet	FI	64
Instruções de serviço	PT	65
Wskazówki bezpieczeństwa	PL	67
Bezpečnostní pokyny	CS	68
Biztonsági utasítások	HU	70
Varnostna opozorila	SL	71
Bezpečnostné pokyny	SK	73
Ohutusjuhised	ET	74
Drošības norādes	LV	76
Nurodymai dėl saugumo	LT	77
Инструкции за безопасност	BG	79
Indicații de siguranță	RO	80
Υποδείξεις ασφάλειας	EL	82

IKA-Werke, Germany  
Reg. No. 004343

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## Contents

	Page
Declaration of conformity	10
Note for USA (FCC)	10
Note for Canada (IC)	11
Operator panel and display	11
Explication of warning symbols	11
Safety instructions	12
Correct use	13
Unpacking	13
Preparations	14
Useful information	15
Commissioning	15
Working with Wireless Controller (WiCo)	15
Interface and output	16
Error codes	16
Maintenance and cleaning	16
Warranty	17
Technical data	17

## Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61326-1, EN 60529 and EN ISO 12100.

Bluetooth® module:

Directive: 1999/5/EC

Standards: EN 300328, EN 301489-1, EN 301489-17, EN 60950-1

## Note for USA (FCC)

This equipment complies with Part 15 of the FCC rules. Any changes or modifications not expressly approved by the Manufacturer could void the user's authority to operate the equipment.

This device complies with Part 15 of the FCC rules subject to the following two conditions:

- (1) This device may not cause harmful interference, and
- (2) This device must accept all interference received, including interference that may cause undesired operation.

### NOTE:

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment.

This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.


## Note for Canada (IC)

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions:

- (1) This device may not cause interference
- (2) This device must accept any interference, including interference that may cause undesired operation of the device.


This device complies with Health Canada's Safety Code 6/IC RSS-210. The installer of this device should ensure that RF radiation is not emitted in excess of the Health Canada's requirement.

## Operator panel and display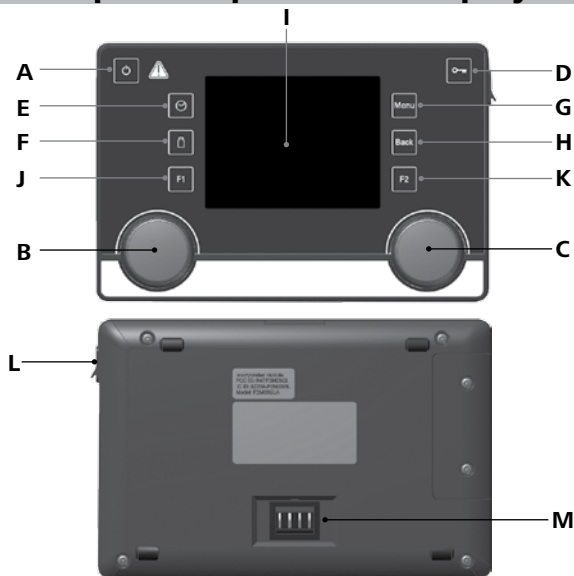

Item	Designation	Function
A	"Power" key:	Turn on/off the WiCo
B	Rotating/pressing knob:	Start/stop the heating function Change the settings of the temperature in working screen
C	Rotating/pressing knob:	Start/stop the stirring function Change the settings of stirring speed in working screen Navigation, selecting and changing the settings in the menu
D	Lock Key:	Lock/unlock knobs and keys
E	Timer button:	Enable timer display
F	Weighing button:	Enable weight display
G	"Menu" key:	Press it once: main menu is displayed Press it a second time: back to the working screen
H	"Back" key:	Return to the previous menu level
I	Display:	Display the settings and actual values
J	F1 key:	Display time/temperature graph
K	F2 key:	Not currently assigned
L	USB interface:	Communication and recharge of WiCo
M	Contacts:	Not currently assigned

## Explication of warning symbols

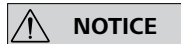Indicates practices which, if not avoided, can result in equipment damage.

# Safety instructions

## Battery pack RB 1

### NOTICE

If during operation the **battery pack RB 1** (rechargeable battery) becomes fully discharged, the device (**Station**) will continue to run or be shut down depending on the value settings for **“Time Out”**, **“Safe Speed”** and **“Safe Temperature”**. If the device is set so that it continues to run when the battery of the **Wireless Controller (WiCo)** is fully discharged, the only means of switching the station off are the **“safe STOP”**, **“ON/OFF”** key and the **“Mains switch”**!

### NOTICE

***Please note the following safety instructions for the battery pack RB 1 (rechargeable battery):***

- Keep the battery pack out of reach of children at all times.
- Store the battery pack in a cool, dry place.
- Never throw the battery pack into a fire. Keep it away from direct sunlight and temperatures above 60 °C. High temperatures will damage the battery pack and render it unusable. Temperatures above 100 °C may cause it to explode.
- Never throw the battery pack into water or expose it to moisture. Water may lead to a short-circuit, causing the battery pack to explode.
- Do not deform or crush the battery pack or damage it in any other way. This can cause battery fluid to leak and/or the battery pack to explode.
- When not in use, keep battery packs away from paperclips,

coins, keys, nails, screws or other small metal objects which could cause the contacts to be bridged. Short-circuiting may result in an explosion.

- Explosion of a battery pack may release battery fluid and cause a fire.
- The lithium polymer battery pack must only be used and charged in **IKA®** products designed for use with this battery pack.
- When the battery pack is inserted it should slide in easily and without resistance. Do not force it.
- If the battery pack is removed for an extended period of time, store it in a sealed plastic bag to prevent short-circuiting due to moisture or contact with metal components.
- The operating temperature range of the battery pack is from 0 °C to + 45 °C. Note that the battery pack capacity will be reduced at temperatures below 20 °C.
- Only the rechargeable battery types recommended in the technical data may be used in the device!

Do not charge batteries that have leaked or that are discolored, deformed or damaged in any other way.

## **Disposal instructions**

- When disposing of the **IKA®** battery pack, please tape over the contacts with adhesive tape to prevent short-circuiting due to moisture or contact with metal components. Short-circuiting may result in an explosion.
- Do not throw used battery packs into your household waste. Dispose of them properly in accordance with statutory regulations.

End users are obliged by law to return all used disposable and rechargeable batteries. Throwing them into the household waste is prohibited. Disposable/rechargeable batteries containing harmful substances are marked with this symbol to indicate that they may not be disposed of as household waste.

- You can return used disposable and rechargeable batteries free of charge to your local authority collection site or to any battery retailer. In doing so you will be complying with statutory regulations and helping to protect the environment.
- Batteries must be disposed of in accordance with local and national regulations.

## Correct use

- **Use**
  - The **Wireless Controller (WiCo)** is used for wireless remote controlling of LR 1000 control and RET control-visc.
- **Range of use (indoor use only)**
  - Laboratories                      - Schools
  - Pharmacies                         - Universities
- **Wireless remote control:**

Before using the wireless link between the **Wireless Controller (WiCo)** and the laboratory device, first check whether your region is included in the radio communication approval for the device. If it is not, remote control can also be performed using a USB cable.

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the device is operated with accessories that are not supplied or recommended by the **IKA®**.
- If the device is operated improperly or in contrary to the **IKA®** specifications.
- If the device or the printed circuit board are modified by third parties.

## Unpacking


- **Unpacking**
  - Please unpack the device carefully;
  - In the case of any damage a fact report must be sent immediately (post, rail or forwarder).
- **Delivery scope**
  - **Wireless Controller (WiCo)**
  - WiCo holder WH 10
  - USB cable type Micro-B/Micro-A (WiCo ↔ device)
  - USB cable type Micro-B/A (WiCo ↔ power supply adapter)
  - Power supply adapter (USB)
  - Screwdriver (Battery cap)
  - Operating instructions
  - A warranty card

## Preparations

- **Charging the Battery Pack RB 1 (rechargeable battery):**


The battery pack in the **WiCo** can be charged by the power supply adapter (included in the delivery).

- **Changing battery to WiCo:**

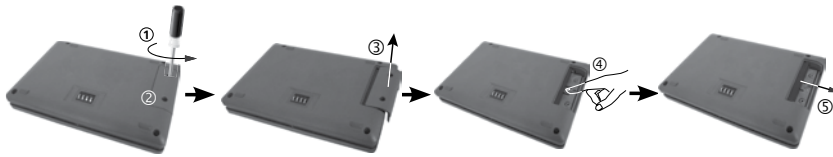Please comply with the relevant safety instructions in the “Safety instructions” section for the RB 1 battery pack!

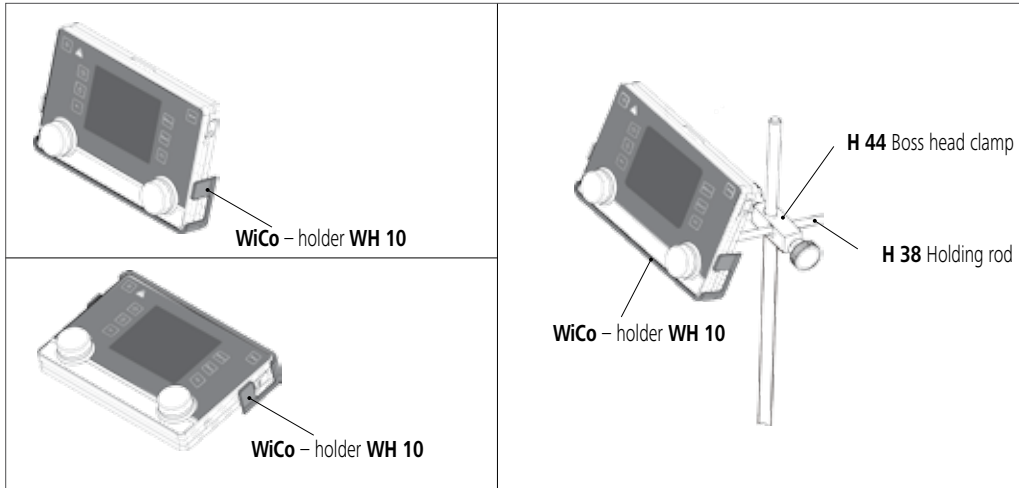- **WiCo – holder WH 10:**

## Useful information


The **WiCo** is equipped with a USB socket (Universal Serial Bus) with which the **WiCo** can be connected with **station**, the USB icon  appears on the screen.

If the **WiCo** is not connected with **station** via a USB cable, the data exchange between the **station** and the **WiCo** via Bluetooth®. In this case, the Bluetooth® icon  is displayed.

Depending on the structure of the building, the **WiCo** can be operated at a distance up to 15 m from the **station**, using the Bluetooth® connection.

## Commissioning

Check whether the voltage given on the type plate corresponds to the available mains voltage.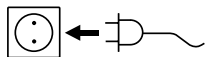

 The socket used must be earthed (fitted with earth contact).

If these conditions have been met, the machine is ready for operation when the mains plug is plugged in.

If these conditions are not met, safe operation is not guaranteed and the machine could be damaged.


Observe the ambient conditions (temperature, humidity, etc.) listed under “**Technical Data**”.

The battery in the **Wireless Controller (WiCo)** must be charged by the power supply adapter before the first use of the device.

The battery could also be charged via the USB port on the **WiCo** (See “**Charging the Battery Pack RB 1**” in “**Preparations**” section).

## First use information (e.g. RET control-visc)

1) Connect **WiCo** to station via USB cable

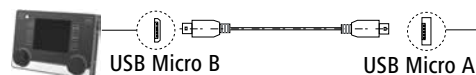- 2) Switch on station and **WiCo**
- 3) Follow the instructions on the **WiCo** screen
- 4) Devices will connect automatically after few seconds
- 5) Remove the USB cable after successful connection for wireless control

## Working with Wireless Controller

After switching on the **WiCo** the operating control changes over from the station screen to the **WiCo** screen. If the **WiCo** is switched off the operating control changes back to the station screen.

### NOTICE

- Emergency-Stop is always possible at the station.
- When operating, please refer to the operating instructions of LR 1000 control and RET control-visc.

## Interface and output

Please refer to the operating instructions of LR 1000 control and RET control-visc.


## Error codes

Please refer to the operating instructions of LR 1000 control and RET control-visc.

## Maintenance and cleaning

The device should operate without the need for maintenance, however it is subject to natural wear and tear on parts and their statistical failure rate.

### Cleaning

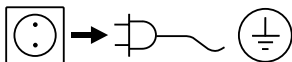Remove the device from the mains before cleaning.

Only clean **IKA®** appliances using these **IKA®** approved cleaning agents.

Dirt	Cleaning agent
Dyes	Isopropyl alcohol
Building materials	Water containing detergent/isopropyl alcohol
Cosmetics	Water containing detergent/isopropyl alcohol
Food	Water containing detergent
Fuels	Water containing detergent
Other materials	Please consult <b>IKA®</b>

- Wear protective gloves during cleaning the instrument.
- Electrical devices may not be placed in the cleaning agent for the purpose of cleaning.
- Do not allow moisture to get into the equipment when cleaning.
- Please consult **IKA®** before using any cleaning or decontamination methods, other than those recommended here.

### Ordering spare parts

When ordering spare parts, please make sure to indicate the following:

- device type
- device serial number; see rating plate
- software version
- position number and description of spare part; see **www.ika.com**.

## Repairs

**Please send in instrument for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.**

For repair, please request the “**Decontamination Certificate**” from **IKA®**, or download printout of it from the **IKA®** website **www.ika.com**.

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

## Technical data

### Wireless Controller

Permitted on time	%	100
Max. communication distance (dependent on the building)	m	15
Dimensions (W x D x H)	mm	160 x 37 x 105
Weight	kg	0.3
Ambient temperature	°C	+ 5 ... + 40
Ambient humidity (relative)	%	80
Protection class according to EN 60529		IP 40
Interface		USB

### RB 1 Battery pack

Voltage	V	3.7
Charging capacity	mAh	2000
Charging time	h	4.5
Working time	h	15
Battery type		Lithium-polymer

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**IKA® - Werke GmbH & Co.KG**