# IKA

designed for scientists

## TW.VX

# IKA

designed for scientists

### IKA-Werke GmbH & Co. KG

Janke & Kunkel-Straße 10, 79219 Staufen, Germany  
Phone: +49 7633 831-0  
eMail: sales@ika.de

#### USA

IKA Works, Inc.  
Phone: +1 910 452-7059  
eMail: sales@ika.net

#### MALAYSIA

IKA Works (Asia) Sdn Bhd  
Phone: +60 3 6099-5666  
eMail: sales.lab@ika.my

#### JAPAN

IKA Japan K.K.  
Phone: +81 6 6730 6781  
eMail: info\_japan@ika.ne.jp

#### VIETNAM

IKA Vietnam Company Limited  
Phone: +84 28 38202142  
eMail: sales.lab-vietnam@ika.com

#### KOREA

IKA Korea Ltd.  
Phone: +82 2 2136 6800  
eMail: sales-lab@ika.kr

#### CHINA

IKA Works Guangzhou  
Phone: +86 20 8222 6771  
eMail: info@ika.cn

#### INDIA

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Discover and order the fascinating products of IKA online: [www.ika.com](http://www.ika.com)

## Explication of warning symbols

### Warning symbols

-  **Danger!** Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.
-  **Warning!** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
-  **Caution!** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
-  **Notice!** Indicates practices which, if not avoided, can result in equipment damage.

### General Symbols

-  Position number
-  Indicates device components relevant to actions.
-  Wrong
-  Indicates the incorrect execution of an action step.
-  Correct / result
-  Indicates the correct execution or the result of an action step.
-  Note
-  Indicates steps of actions that require particular attention.

## Safety instructions

### General information

- > Read the operating instructions in its entirety before using the device and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.
- > Keep the operating instructions in a place where it can be accessed by everyone.
- > Ensure that only trained staff work with the device.
- > Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

### Device design

-  **Notice!**
- > Keep the adhesive surface free of dust.
- > Check the adhesive force of the adhesive pad after each removal and replacement.
- > Clean the adhesive pad of the shaking attachment with a surfactant-containing water moistened, lint-free cloth.
- > Protect the device and accessories from bumps and impacts.
- > Check the device and accessories for damage before each use. Do not use damaged components.

### Working with the device

-  **Danger!**
- > To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
-  **Warning!**
- > Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- > Beware of hazards due to:
  - flammable materials,
  - glass breakage as a result of mechanical shaking power,
  - incorrect container size,
  - overfilling of media,
  - unsafe condition of container.
- > Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
  - Keep the sample container as vertical as possible.
  - Only exert the minimum pressure necessary for the mixing process.
  - Use attachments for multiple samples when there is a large number of samples.
- > If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
- > Reduce the speed if:
  - the medium splashes out of the vessel because the speed is too high,
  - the appliance is not running smoothly,
  - the appliance begins to move around because of dynamic forces.
- > Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.
- > Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
- > Sharp edged vessels cause wear debris on the attachments.
- > When handling hazardous substances or mixtures of substances which are toxic or are contaminated with pathogenic micro-organisms, the user should take appropriate precautions. Observe the national regulations, the biosafety level of your laboratory, the material safety data sheets.
- > For germs or biological material or risk group II or higher, please observe the "Laboratory Biosafety Manual" issued by the WorldHealth Organization).
-  **Caution!**
- > Wear your personal protective equipment in accordance with the hazard category of the media to be processed. There may be a risk from:
  - splashing and evaporation of liquids,
  - ejection of parts,
  - release of toxic or combustible gases.

### Accessories

- > Safe operation is guaranteed only with the use of original IKA accessories.
- > Observe the operating instructions of the accessories.

### Disposal instructions

- > The device, accessories and packaging must be disposed of in accordance with local and national regulations.

## Intended use

### Use

- > The TW.VX is a shaking attachment for mixing liquids.
  - Touch mode for single test tubes.
  - Continuous mode for single or several vessels.

### Area of use

- > Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- > The safety of the user cannot be guaranteed:
  - if the device is operated with accessories that are not supplied or recommended by the manufacturer,
  - if the device is operated improperly or contrary to the manufacturer's specifications,
  - if the device or the printed circuit board are modified by third parties.

## Unpacking

### Unpacking

- > Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

### Scope of delivery

1	TW.VX	4	MS 1.21 One-hand insert
2	MS 3.1 Standard attachment	5	Lever tool
3	MS 3.3 Universal attachment	6	Operating instructions

## Assembly

1. Degrease the installation surface of the drive unit (e.g. with isopropanol).
2. Remove the protective film from the adhesive mat.
3. Place the shaking attachment in the center of the TWISTER drive unit.
4. Remove the shaking attachment with the lever tool, by lifting it upwards.

## Cleaning

- > Use only cleaning agents which have been approved by IKA to clean IKA devices. These are: water (containing surfactant) and isopropyl alcohol.
- > Wear protective gloves during cleaning the devices.
- > Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.
- > The adhesive force of the adhesive pad can be restored by cleaning it with water.

## Accessories

- > Permitted speed ranges of the attachments:

Speed range (rpm)		
0/100 ... 2500	0/100 ... 1300	0/100 ... 500
MS 3.1	MS 3.3, MS 3.5, MS 3.51, MS 1.21, MS 1.30, MS 1.31, MS 1.32, MS 1.33, MS 1.34, MS 3.4	STICKMAX II

- > For accessories see [www.ika.com](http://www.ika.com).

## Technical data

Type of movement	orbital	Dimensions (W x H)	94 x 28 mm
Shaking stroke	4.5 mm	Weight	0.54 kg
Permissible shaking weight (incl. attachment)	0.5 kg		

> Subject to technical changes!