designed for scientists

Montageanleitung / Assembly instructions I-MAG cart

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Read the assembly instructions in its entirety before using the I-MAG cart and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.

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Safety instructions

Caution!

- › Ensure that your fingers do not get crushed when setting down the I-MAG cart.

Caution! **Notice!**

- › Set up the I-MAG cart in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- › Observe the maximum permissible load-bearing capacity. (see “Technical data”)
- › Beware of the high dead weight of the I-MAG cart when transporting.
- › The center of gravity of the vessel must lie in the center of the I-MAG cart.
- › Wear protective gloves during assembling the I-MAG cart.
- › Do not use your hands to lock / unlock the castors.
- › Do not start stirring if the castors are unlocked or the cart is moving.
- › Do not start stirring if the I-MAG cart is on a ramp or an uneven floor.

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Unpacking

1	Chassis assembly
2	Handle bar
3	Holding plate
4	M10 x 25 hexagon socket screw and screw washer (2x)
5	M5 x 35 hexagon socket head screw (2x)
6	Cable clip (5x)
7	ISO 2936-4 Screwdriver
8	ISO 2936-8 Screwdriver

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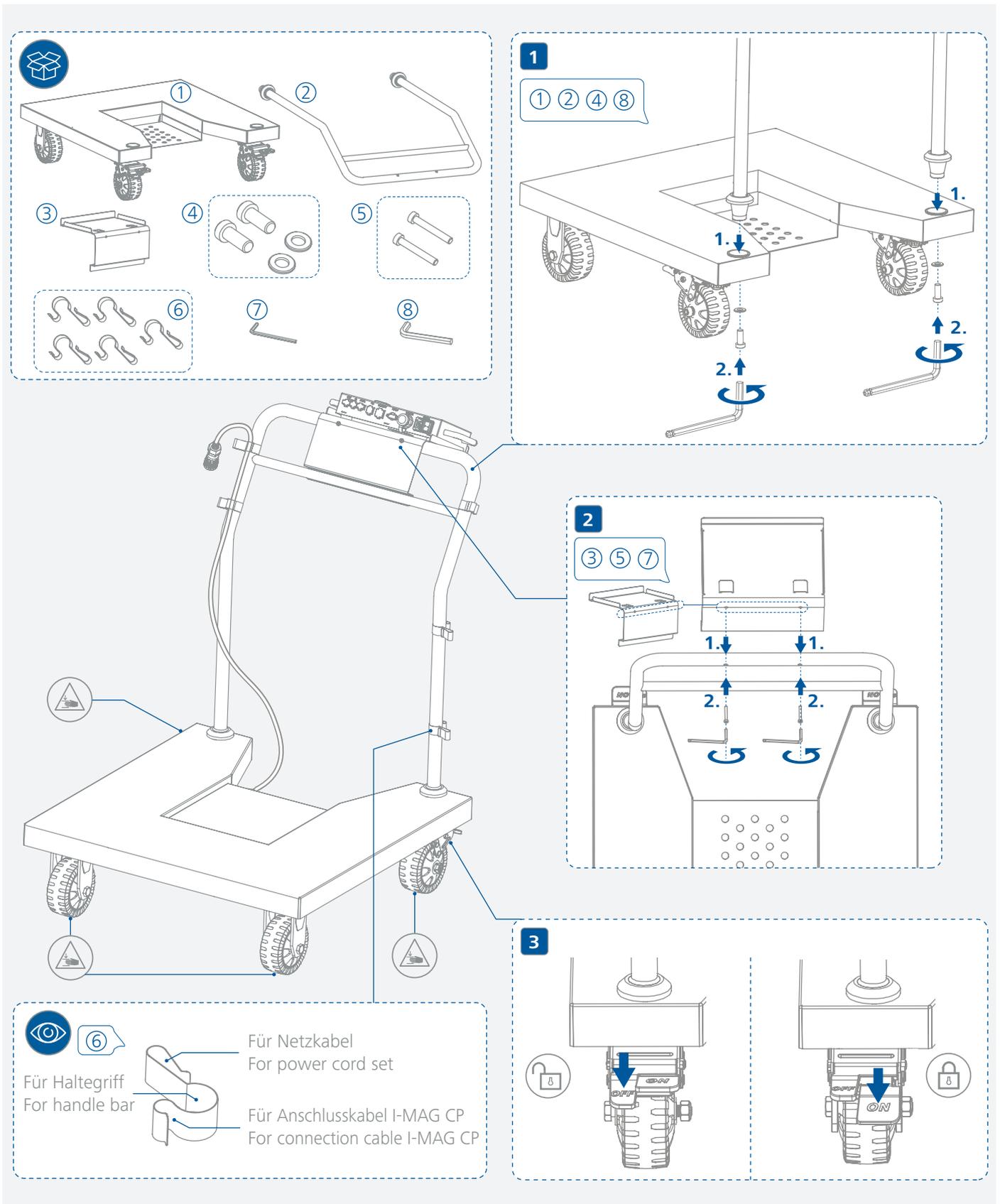
Technical data

Dimensions (W x D x H)	700 x 995 x 970 mm
Weight	32 kg
Bearing load Max.	400 kg
Seating plate material	Stainless steel 1.4301
Castor tread material	PA

Subject to technical changes!

IKA-Werke GmbH & Co. KG

Janke & Kunkel-Straße 10, 79219 Staufen, Germany • Phone: +49 7633 831-0 • Fax: +49 7633 831-98 • eMail: sales@ika.de

Discover and order the fascinating products of IKA online: www.ika.com

Technical specifications may be changed without prior notice.

20000037826_20116007_I-MAG cart_052022