
IKA

designed for scientists

C 6010/C 6012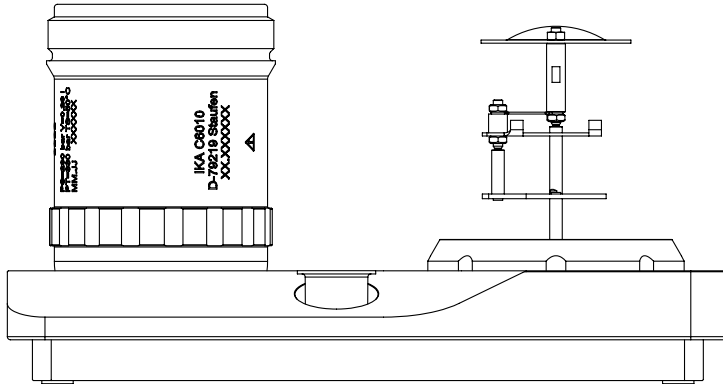

Betriebsanleitung	DE	2
Ursprungssprache		
Operating instructions	EN	11
Mode d'emploi	FR	20
Руководство пользователя	RU	29
使用说明	ZH	38
Indicaciones de seguridad	ES	47
Veiligheidsinstructies	NL	49
Norme di sicurezza	IT	51
Säkerhetsanvisningar	SV	53
Sikkerhedshenvisninger	DA	54
Sikkerhetsinformasjon	NO	56
Turvallisuusohjeet	FI	58
Instruções de segurança	PT	59
Wskazówki bezpieczeństwa	PL	61
Bezpečnostní upozornění	CS	63
Biztonsági utasítások	HU	65
Varnostna navodila	SL	67
Bezpečnostné pokyny	SK	68
Ohutusjuhised	ET	70
Drošības norādes	LV	72
Saugos nurodymai	LT	74
Инструкции за безопасност	BG	75
Indicații de siguranță	RO	77
Υποδείξεις ασφαλείας	EL	79

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Source language: German

Contents

	Page
Declaration of Conformity	11
Explanation of symbols	11
Safety instructions	11
Included with delivery C 6010/C 6012	13
Opening the decomposition vessel	14
Fastening cotton thread	14
Closing the decomposition vessel	15
Combustion procedure	15
Cleaning the decomposition vessel	15
Halogens C 6012	15
Care and maintenance	16
Replacing the ignition wire	16
Replacing the ignition electrode	17
Replacing the crucible holder	17
Replacing the ground electrode	17
Maintaining Oxygen valve	17
Replacing C 6010 sealing ring	18
Leakage test	18
Degas	18
Decomposition vessel identification	19
Technical data	19
Warranty	19
Spare parts list	94
Spare parts diagram	95

Declaration of Conformity

We declare under our sole responsibility that the product to which this declaration relates is in conformity with the Pressure Vessel Directive 2014/68/EU and the regulations AD 2000. A copy of the complete EU Declaration of Conformity or further declarations of conformity can be requested at sales@ika.com.

Explanation of symbols

This symbol identifies information **that is absolutely essential for your safety and health.** Failure to observe this information may result in negative effects for your health and injuries.

This symbol identifies information that is important for problem-free operation as well as for working with the decomposition vessel. Failure to observe these instructions may result in inaccurate measurement results or damage to the decomposition vessel.

Safety instructions

Intended purpose:

The C 6010 and C 6012 decomposition vessels must only be used for determining the gross calorific value of solid and liquid combustible materials in accordance with DIN 51900, BS 1016 T5, ISO 1928, ASTM 5468, ASTM 5865, and ASTM 4809. Operation of decomposition vessels is only permitted in combination with the IKA calorimeter C 6000 g/iso.

Operating requirements:

The maximum **amount of energy** input into the decomposition vessel must not exceed **40000 J**. (Select the weight of the sample accordingly). The permissible **operating pressure of 230 bar (23 MPa)** must not be exceeded. The maximum permissible **operating temperature** must not exceed **50 °C**. Do not fill the decomposition vessel too full of the sample. Only fill the decomposition vessel with oxygen up to a maximum pressure of **40 bar (4 MPa)**. Monitor the adjusted pressure on the pressure reducer.

Perform a check before every combustion to ensure there are no leaks (see sections "**Leakage test**").

Explosive substances:

Many substances tend to combust in an explosive manner (for example because of the formation of peroxide). This may cause the decomposition vessel to burst.

The C 6010 and C 6012 standard decomposition vessels must not be used for examinations on samples that are capable of exploding. It is absolutely essential to use a special high-pressure decomposition vessel to contain the sample in these cases!

Notes on the sample:

Substances of which the combustion behavior is not known must be examined for their combustion behavior before combustion in the decomposition vessel C 6010 and C 6012 (danger of explosion).

If you are burning **unknown samples**, leave the room or **keep a safe distance** between you and the calorimeter.

Benzoic acid must only be burned in the form of pellets! Combustible dust and powder must be compressed into pellets before combustion. Oven-dry dust and powder such as wood chips, hay, straw, etc. burn in an explosive manner! They must be moistened first! Readily combustible liquids with a low vapor pressure must not be come in direct contact with the cotton thread (for example tetramethyl dihydrogen disiloxan)!

Combustion residue, auxiliary materials:

In addition, toxic residues of combustion are possible in the form of gasses, ash or precipitates on the inner wall of the decomposition vessel, for example.

Observe the accident prevention requirements applicable to the activity and the work station. Wear personal safety equipment.

When handling combustion samples, combustion residues and auxiliary materials, the appropriate safety requirements must be observed. The following are examples of substances that may cause dangers:

- corrosive
- easily flammable
- capable of exploding
- contaminated with bacteria
- toxic

Oxygen:

When working with oxygen, observe the appropriate requirements.

Danger warning: As a compressed gas, oxygen promotes combustion, supports combustion intensively and may react violently with combustible substances. **Do not use any oil or grease!**

Using a crucible made of stainless steel:

When using crucibles made of stainless steel, their condition should be carefully checked after every experiment.

A reduction in the thickness of the material may cause the crucible to burn and may damage the C 6010 and C 6012 decomposition vessel.

For reasons of safety, crucibles must not be used any more after a maximum of 25 combustion procedures.

Ignition wire:

The decomposition vessel C 6010 and C 6012 are equipped with a fixed ignition wire. Since the ignition wire is subjected to a high degree of material stress during combustion, it should be replaced after 100 tests for C 6010 (kanthal) and after 50 tests for C 6012 (platinum) at the latest.

Specification of the decomposition vessel:

The decomposition vessel is manufactured in accordance with the regulation for pressure vessels **2014/68/EU and the regulations AD 2000**. This can be recognized from the **CE symbol** with the identifying number of the testing station named. The decomposition vessel is a pressure device of Category III. The decomposition vessel has been subjected to an EC prototype test. The CE declaration of conformity represents our guarantee to you that this decomposition vessel complies with the pressure device described in the EC prototype test certificate. The decomposition vessel has been subjected to a pressure test at a **test pressure of 330 bar (33 MPa)** and a leak test with oxygen at 30 bar (3 MPa).

C 6010 and C 6012 decomposition vessels are **experiment autoclaves** and must be tested by a professionally trained person each time before they are used.

An individual application is understood here to mean a series of experiments that are performed under roughly the same conditions in terms of pressure and temperature. Experiment autoclaves must be operated in special chambers (C 6000 gs/iso).

Repeated tests:

The decomposition vessel must be subject to repeated tests (internal tests and pressure tests) by a **person with professional training**. The intervals between tests must be determined by the operator based on experience, operating manner and the material used in the decomposition vessel.

The declaration of conformity loses its validity if mechanical modifications are made to the experiment autoclaves or if stability can no longer be guaranteed as a result of heavy corrosion (e.g. pitting by halogens, see section "Halogens C 6012").

The **threading** on the body of the decomposition vessel and cap screw in particular are subject to a high level of mechanical stress and must therefore be monitored regularly for **wear and tear**.

The condition of the seals must be checked for functionality must be ensured by means of a test for leaks (see sections "**Leakage test**").

Pressure tests and service tasks on the decomposition vessel must only be performed by **persons with professional training**.

We recommend that the decomposition vessel be sent into our factory for inspection and repairs if necessary after either 1000 experiments or after one year or, depending on the application, even sooner than this.

Definition of person with professional training:

A person with professional training as defined in these operating instructions is someone:

1. whose training, knowledge and experience gained through practical activities ensures that that person will perform the tests in a proper manner.

2. who is sufficiently reliable.
3. who is not subject to any instructions in terms of testing activity.
4. who is equipped with suitable testing equipment if necessary.
5. who can provide suitable proof demonstrating compliance with the requirements listed in 1.

Operating pressure containers:

National regulations and laws for operating pressure containers must be observed! Anyone who operates a pressure container must maintain it in proper condition, must monitor it and perform necessary maintenance and repair tasks without delay, and must take measures appropriate for the circumstances to ensure safety.

A pressure container must not be operated if it exhibits defects that could endanger those working with it or third parties.


You can obtain a copy of the pressure vessel regulation from Beuth Press.

Use only original spare parts from IKA!


Included with delivery C 6010/C 6012

- 1 x Decomposition vessel
- 1 x Quartz dish C 4
- 1 x Operating instructions
- 1 x Emergency-vent knob
- 1 x SW 8/10 double open ended wrench
- 2 x SW 5,5/7 double open ended wrench
- 1 x SW 17 Wrench
- 5 x Kanthal ignition wire (C 6010)
- 2 x Platinum ignition wire (C 6012)
- 1 x O-ring set
- 1 x C 710.4 cotton thread
- 1 x Test report

Opening the decomposition vessel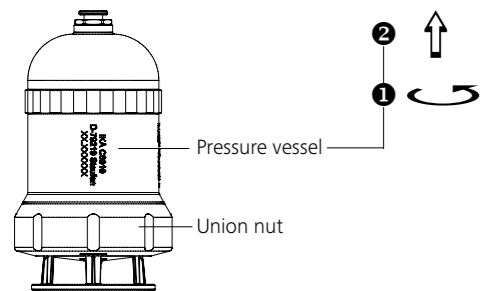


The decomposition vessel consists of the following individual parts: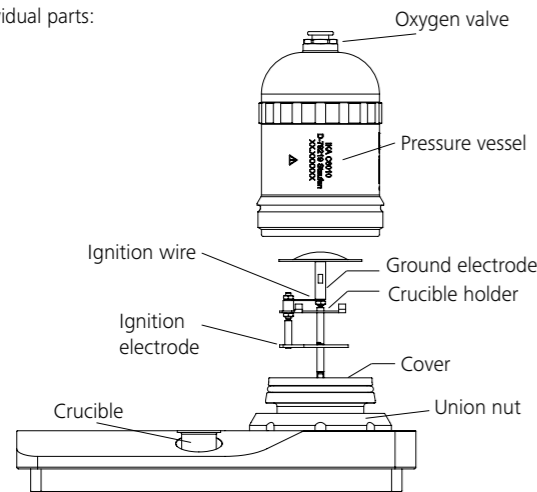


Fastening cotton thread

Note:

The straightening and assembly work is done with the help of C 6010 organizer.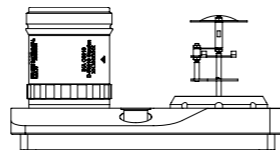


The decomposition vessel is equipped with a fixed ignition wire. A cotton thread has to be fastened to the ignition wire in order to carry out combustion.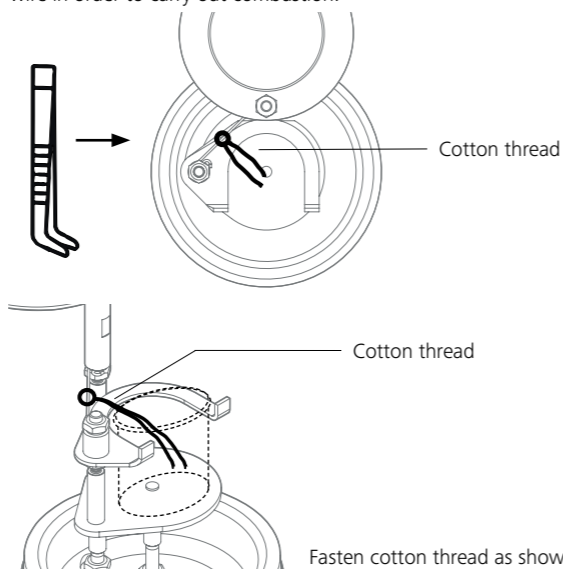

Fasten cotton thread as shown in the images.

Closing the decomposition vessel

- Place the cover into the union nut.
- Screw the pressure vessel into the union nut until stop.

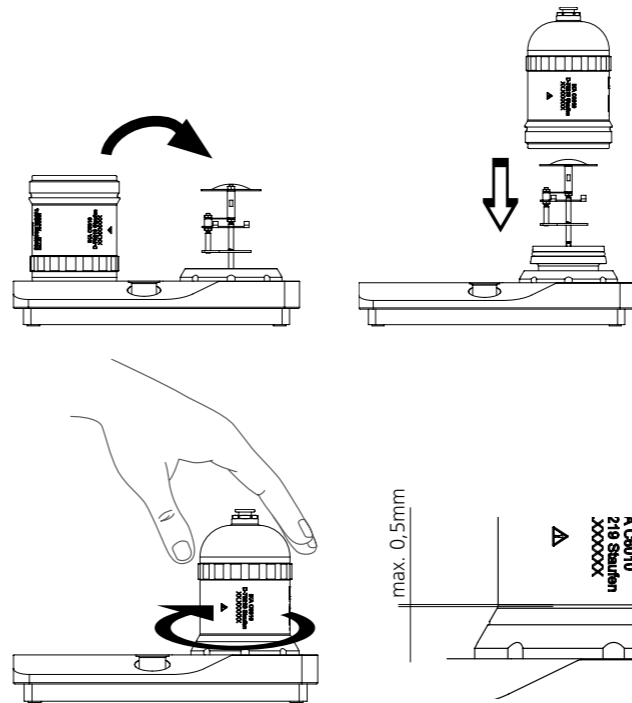Combustion procedure

For the decomposition vessel charging procedure with fuel samples and the general procedure for performing combustion please refer to the relevant standards and the operating instructions for the IKA C 6000 g/iso. When working with C 6000, the oxygen filling and degassing to the decomposition vessel is done inside the device.

Cleaning the decomposition vessel

After each combustion the condensate has to be cleaned from all of the parts of the decomposition vessel which come into contact with the combustion products. It suffices to wipe out the parts with an absorbent, non-fibrous cloth. Combustion residues in the crucible, e.g. carbon black or ash are removed in the same manner. The decomposition vessel may not be mechanically cleaned under any circumstances if it is not capable of being cleaned in the manner described (e.g. due to burn-in or pitting).

In such cases please contact your **competent service center** or send in the decomposition vessel to the factory for inspection and cleaning.

Halogens C 6012

Pitting which severely impairs operator safety may be caused on the inner surfaces of the decomposition vessel from combustion products which contain halogen. If the halogen content of the samples is significantly greater than the trace range, then decomposition vessel C 6012, which is specially designed for such applications, shall be used.

To enable the **quantitative evaluation of halogens and sulphur**, the C 6012 decomposition vessel is preconditioned before shipment. This means that the inner surfaces and the fittings of the decomposition vessel are catalytically activated by means of a patented procedure.

Parts thus treated are recognizable because of their opaque, tarnished and partly spotty surfaces. This appearance of this surface is intentional and should not be confused with traces of wear.

This state is important for the following analysis and may not be treated with brushes or other hard materials.

Wiping and drying should be done using a soft cloth.

The C 6010 decomposition vessel can not be used for such applications because of the safety risk and insufficient analysis results.

Care and maintenance

The work described in the following may only be performed by personnel with professional training or **IKA**.

All work on the fittings of the decomposition vessel require a leak test and new pressure testing.

Furthermore, reference is made to the regulation for pressure Vessels 2014/68/EU.

Send in the case of service the equipment back in the original packaging. Storage packing is not sufficient for the back dispatch. Use additionally a suitable packing.

Replacing the ignition wire

The C 6010 and C 6012 decomposition vessels are equipped with a fixed ignition wire. Since the ignition wire is subjected to a high degree of material stress during combustion, it should be replaced after 100 tests for C 6010 (kanthal) and after 50 tests for C 6012 (platinum) at the latest. However, this largely depends on the type of samples fired.

The ignition wire must also be replaced if corrosion occurs.

NOTE

The crucible holder must not be twisted while doing so. Leakages of the decomposition vessel are the results (see the section "Leakage test").

Leakages at the decomposition vessel container can cause heavy damage to the calorimeter and the decomposition vessel. Ignoring can lead to health impairment and injury.



Insert/Tightening of the ignition wire:

The ignition wire is shaped in such a way that the electrodes of the decomposition vessel fit into the eyes.

The following points have to be strictly observed:

- Align the points horizontally between the electrodes
- Do not press the electrodes together. The shielding of the ignition electrodes (13) and the crucible holder (item 8) must not touch the ground electrode (item 7). (see details "Z")
- The ignition electrode must not be twisted and thereby loosened
- Follow the sequence in reverse to tighten the nut
- **Ensure that there is a sound contact with the ignition wire.**


Loosen the ignition wire (item 41) from ground electrode (item 7) and crucible holder (item 8).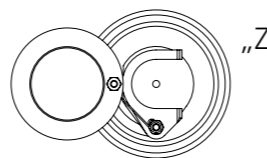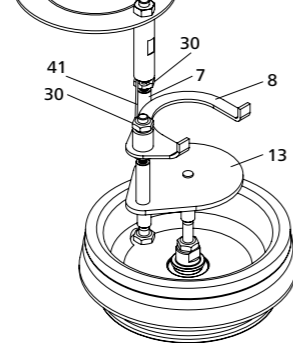

- Item 7 Ground electrode
- Item 8 Crucible holder
- Item 13 Ignition electrode
- Item 30 Hex nuts M3
- Item 41 Ignition wire


Replacing the ignition electrode

Take care two new o-rings (item 10 and 23) when inserting the new electrode.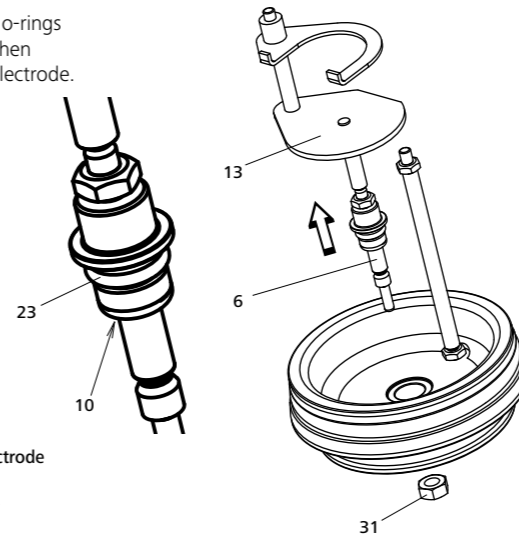

- Item 6 Shaft
- Item 10 O-ring
- Item 13 Ignition electrode
- Item 23 O-ring
- Item 31 Hex nut M4

Replacing the crucible holder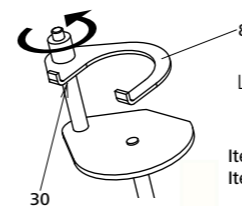


Loosen hex nut (item 30).

- Item 8 Crucible holder
- Item 30 Hex nut M3

Replacing the ground electrode

Loosen hex nut (item 30).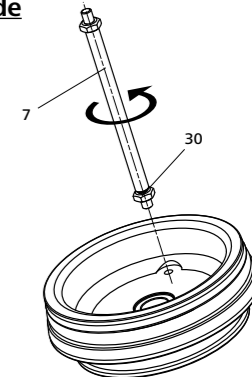


- Item 7 Ground electrode
- Item 30 Hex nut M3

Maintaining oxygen valve


The oxygen valve is used to fill and empty the decomposition vessel. The o-ring (item 22) may wear because of aggressive combustion gases, which means that the decomposition vessel may not be tight.

The oxygen valve must be unscrewed to change the o-ring.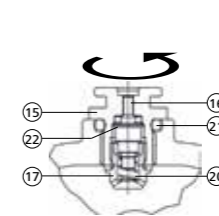


Mounted oxygen valve:

- Item 15 Screwing
- Item 16 Valve piston
- Item 17 Nozzle
- Item 20 Spring
- Item 22 O-ring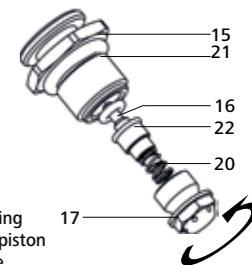


Replacing C 6010 sealing ring

Remove the seal ring (item 11) as shown in the following images.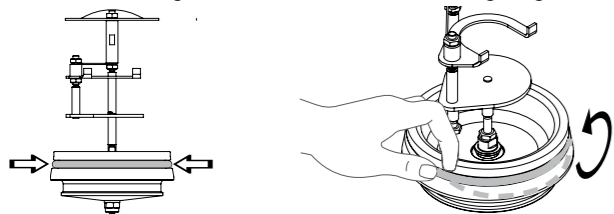

The new sealing ring has to snap exactly as indicated in following images.

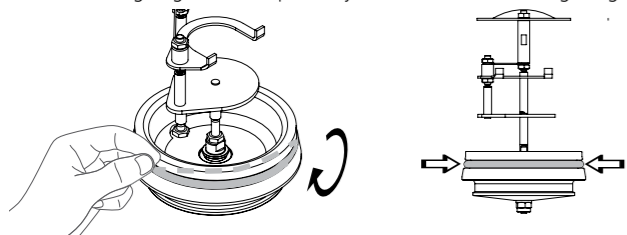A leak test and pressure testing shall be performed after assembly has been successfully completed (see the section "Leakage test").

Leakage test

Perform a leakage test before every combustion process. Always check the tightness of decomposition vessel after carrying out maintenance.


Follow the safety instructions.

- Fill the decomposition chamber in the calorimeter system or using the C 248 oxygen filling station
- Remove the decomposition vessel and immerse fully in a container filled with water
- Check the decomposition vessel for oxygen leaks by watching for bubbles
- If bubbles are seen then the decomposition vessel is not sealed correctly.
- Vent the decomposition vessel and check the seals

Never use a leaking decomposition vessel in the calorimeter.

Degas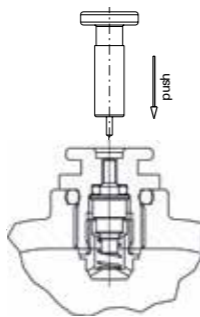


For the C 6000, venting of the decomposition vessel takes place inside the device. The vent knob is supplied for manual release of tension in the decomposition vessel. (Even in case of error function of the filling unit in the calorimeter C 6000 iso / C 6000 gs)

Escaping decomposition gases might be detrimental to health. Please work under a fume hood.

Decomposition vessel identification

The individual parts of the decomposition vessel are clearly identified by means of the engraved serial number and the technical data.

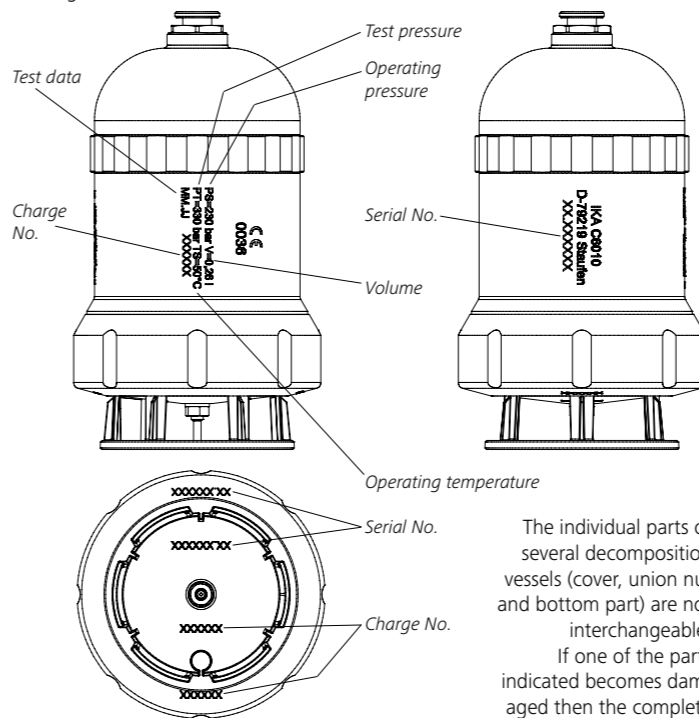The individual parts of several decomposition vessels (cover, union nut and bottom part) are not interchangeable.

If one of the parts indicated becomes damaged then the complete decomposition vessel has to be sent to the factory for testing.

Technical data

Operating pressure	PS	max. 230 bar
Test pressure	PT	330 bar
Volume	V	260 ml
Operating temperature	TS	max. 50 °C
Sample decomposition		isochore
Oxygen valve		plate valve
(vacuum operation not possible)		
Crucible holder		as in DIN 51900
Material in contact with the product		
C 6010		1.4462
C 6012		2.4610.40

Subject to technical changes!

Warranty

In accordance with IKA Terms and Conditions of Sale, the warranty period is 12 months. For claims under the warranty please contact your local dealer. You may also send the device direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited
IKA Works (Thailand) Co. Ltd.