IKA

designed for scientists

C 6030

Betriebsanleitung	DE	4
Ursprungssprache		
Operating instructions	EN	6
Mode d'emploi	FR	10
Руководство пользователя	RU	13
使用说明	ZH	17

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DE

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Source language: German

Contents

	Page		Page
EU Declaration of conformity	2	Setting up of the instrument	8
Contents	6	Function	8
Explanations of symbols	6	Maintenance and cleaning	9
Safety instructions	6	Fault table	9
Unpacking	7	Warranty	9

EN

Explanations of symbols

General hazard

This symbol identifies information **which is of absolute importance for the safety of your health**. Non-compliance may lead to impairments to health and injury.

This symbol identifies information which is important for proper operation and handling of the decomposition vessel. Non-compliance may result in imprecise measurements or damage to the venting station.

Safety instructions

Read the operating instructions in full before starting up and follow the safety instructions.

Keep the operating instructions in a place where they can be accessed by everyone.
Ensure that only trained staff work with the appliance.

Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

6

Furthermore, toxic combustion residue in the form of gases, ashes or condensation, for example, is possible in the inner wall of the washing bottle.

Observe the accident prevention requirements applicable to the activity and the work station.

Wear your personal protective equipment.

When handling combustion samples, combustion residue and auxiliary materials, please observe the relevant safety regulations. The following materials, for example, could pose a risk:

- corrosive
- highly flammable
- explosive
- bacteriologically contaminated
- toxic

Combustion gases are hazardous to health, therefore the venting hose must be connected to a suitable gas cleaning system or extraction system.

The C 6030 venting station is used for controlled venting of **IKA** decomposition vessels C 6010 and C 6012 in conjunction with the **IKA** C 6000 calorimeter system in the "decomposition" operating mode.

The venting speed may be infinitely adjusted via a throttle valve by using a manual adjusting screw. Please note that the throttle valve is **closed** at the beginning of the venting operation.

Do not operate without the ground clamp included in the delivery quantity!

The maximum permitted operating pressure is 40 bar (at the vent header **(f)**).

Please observe the general laboratory guidelines when handling glass vessels.


Unpacking

• Unpacking

- Please unpack the device carefully
- In the case of any damage a report must be sent immediately (post, rail or forwarder).

• Delivery quantity


- Base plate
- Gas washing bottle
- GL screw joint
- Pressure tubing
- Drainage hose
- Venting head
- Fork spanner
- Set of O-Rings
- Operating instructions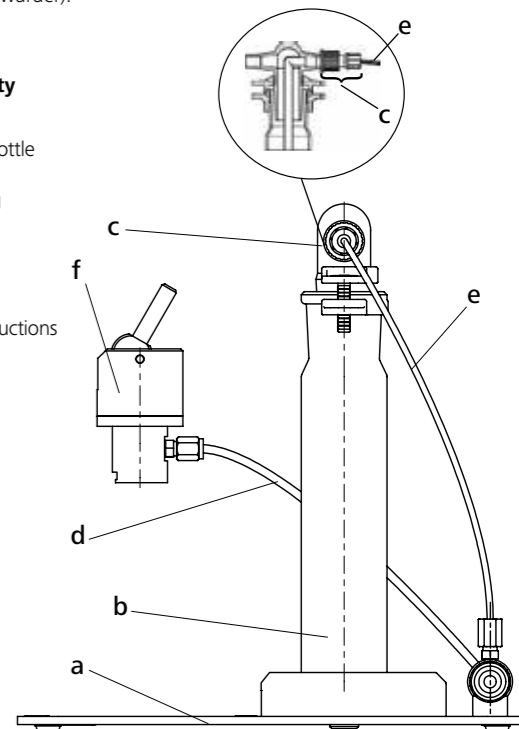

7


Setting up of the instrument

Place the venting station on a horizontal surface and insert the gas washing bottle into the gas washing bottle holder until it comes to a stop. Connect the drainage hose (e) to the throttle valve and the GL screw joint of the washing bottle as shown in the diagram. The tubing is fastened by closing the GL screw joint.

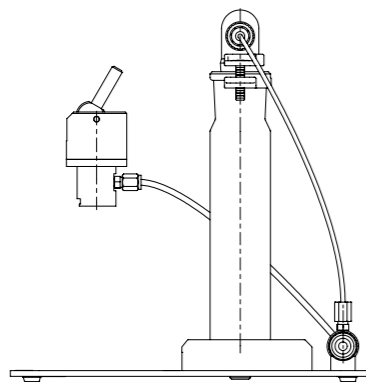WARNING Now fasten the pressure tubing (d) to the venting head and the throttle valve.

All of the screw joints are tightened by means of the A/F 10 fork spanner (1/4 turning) which is included in the delivery quantity.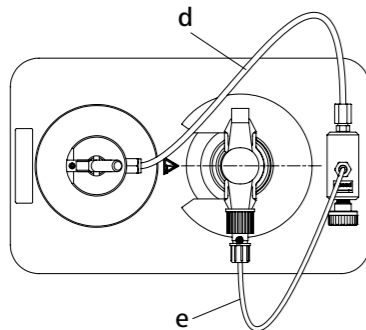

Function

Pour an absorption solution that is suited to your purposes into the gas washing bottle and **secure the ground joints by means of the ground clamp**. Close the throttle valve by slightly turning the adjusting screw in a clockwise direction.

Note: There may be some formation bubbles in the system when the adjusting screw is hand-tightened with maximum operating pressure in the decomposition vessel.

Remove the decomposition vessel from the calorimeter and place it into the venting station. Push the venting head onto the decomposition vessel until it comes to a stop. The valve of the decomposition vessel is opened by pressing down the self-locking eccentric lever.

The venting head cannot be separated from the decomposition vessel in this state.

Check to make certain that the ground clamp fits correctly onto the gas washing bottle and slowly open the throttle valve.

Take note of any bubble formation and correct in order to achieve the venting speed desired.

The venting head is removed when the decomposition vessel is completely expanded (lever in upward position) and the throttle screw is closed again.

WARNING The gas washing bottle may be damaged if venting is performed too quickly.

Maintenance and cleaning

Clean the capillary boring of the venting flask from time to time through expansion with a decomposition vessel filled with pure oxygen-filled without a gas washing bottle attached.

Check the condition of the sealing washer on the filling flask and replace it if necessary.

Flask replacement / removing the pressure spring: The venting head is disassembled by loosening the three countersunk screws located on it. The flask can then be completely removed.

Use only cleaning agents which have been approved by IKA to clean **IKA** devices. These are water (with tenside) and isopropanol.

- Do not allow moisture to get into the appliance when cleaning.
- Wear protective gloves when cleaning the devices.
- Please consult with **IKA** before using any cleaning or decontamination methods, other than those recommended here.

Ordering spare parts

When ordering spare parts, please give:

- Device type
- serial number, see type plate
- Item number and designation of the spare part, see www.ika.com.

Repair

Please send instrument in for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For this you should request the "Decontamination Certificate" from **IKA**, or use the download printout of it from the **IKA** website www.ika.com.

Return the instrument in its original packaging. Storage packaging is not sufficient. Also, please use suitable shipping package materials.

Fault table

Error:	Decomposition valve is not expanded.
Cause:	Flask capillary is deformed.
Remedy:	Replace flask.
Error:	Leakage on venting head when expanding.
Cause:	Sealing washer missing or defective.
Remedy:	Replace sealing.
Error:	Flask in venting head does not retract.
Cause:	Flask boring is dirty or the pressure spring is defective.
Remedy:	Clean or replace spring.
Error:	Leakage on throttle valve
Cause:	O-ring on throttle screw defective.
Remedy:	Replace O-ring.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 12 months. For claims under the warranty please contact your local dealer. You may also send the machine directly to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited