2 x Maulschlüssel DIN 894 SW 12
 2 x Open end wrench DIN 894 A/f 12
 2 x Clé plate DIN 894 ouverture 12

Fig. 2

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Assembling the dispersion tool (overview fig. 3)

- Open the knurled screw (item 3) so that the thread does not protrude into the hole.
- Insert the dispersing element (item 2) into the drive unit (item 1) as far as the stop. After some resistance (depending on the model of the drive - pressure ball) you will hear the shaft clicking into place. The dispersing element is fitted correctly if the recessed edge of the dispersing element is flush with the bottom edge of the drive unit.

- **Attention:** Now secure the dispersing element by screwing the knurled screw tight (item 3).

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IKA Japan K.K.
 Phone: +81 6 6730 6781
 eMail: info_japan@ika.ne.jp

INDIA

IKA India Private Limited
 Phone: +91 80 26253 900
 eMail: info@ika.in

UNITED KINGDOM

IKA England LTD.
 Phone: +44 1865 986 162
 eMail: sales.England@ika.com

VIETNAM

IKA Vietnam Company Limited
 Phone: +84 28 38202142
 eMail: sales.lab-vietnam@ika.com

THAILAND

IKA Works (Thailand) Co. Ltd.
 Phone: +66 2059 4690
 eMail: sales.lab-thailand@ika.com

TURKEY

IKA Turkey A.Ş.
 Phone: +90 216 394 43 43
 eMail: sales.turkey@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com

IKAworldwide

IKAworldwide /// #lookattheblue

@IKAworldwide

Technical specifications may be changed without prior notice.

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Safety instructions					
<ul style="list-style-type: none">Read the operating instructions in full before starting up and follow the safety instructions. Keep the operating instructions in a place where they can be accessed by everyone. Ensure that only trained staff work with the appliance. Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations. Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of splashing and evaporation of liquids; body parts, hair, clothing and jewellery getting caught. No liquid may get into the drive unit of the disperser. The work has to be interrupted immediately if you notice unusual noise. The bearings and seals must be controlled, and replaced if necessary. Note the operating instructions of the disperser unit. Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on. Do not use any damaged dispersion tools e.g. with hairline cracks or splits. <div><div><div>⚠ Risk of injury! When working with the R 50 agitator shaft the safety cage must not be removed. The agitator shaft must be installed in the T 50 drive unit such that the bent safety cage prevents any direct contact with the shaft from in front.</div></div></div> <div><div><div>⚠ The bearing bushing can be very hot during the operation. Let the dispersing tool cool down after the operation before touch and remove it from the device.</div></div></div>					
Product information					
<ul style="list-style-type: none">The R 50 stirring shaft is a high-speed ball-bearing rotary shaft. A splash guard seals it against the medium. The maximum immersion depth of the dispersion tool is 180 mm. The free-wheeling conical shaft is dangerous. Observe the safety instructions! Content of package: R 50 stirring shaft with dissolver disc R 1402 and tool. The generator (dispersion head) consists of the rotor R 1405 propeller made of stainless steel. This part can be unscrewed from the shaft.					
Application instruction					
<ul style="list-style-type: none">Remove the protection cap before use. Use this cap to protect shaft tubes from damage when not in service. Dispersing causes the medium to heat up. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. The function of the dispersion tools depends on the condition of the sharp edges of the rotor. Abrasive media can round off these edges fast, whereby the disperse effect diminishes.					
Designation	Application	Max. permissible speed [rpm]	Working range pressure [bar]	Working range vacuum [mbar]	
R 50	Wet crushing, suspension, emulsifying, deagglomerating (Batch-operation)	10,000	-	-	

Product information					
<ul style="list-style-type: none">The R 50 stirring shaft is a high-speed ball-bearing rotary shaft. A splash guard seals it against the medium. The maximum immersion depth of the dispersion tool is 180 mm. The free-wheeling conical shaft is dangerous. Observe the safety instructions! Content of package: R 50 stirring shaft with dissolver disc R 1402 and tool. The generator (dispersion head) consists of the rotor R 1405 propeller made of stainless steel. This part can be unscrewed from the shaft.					
Application instruction					
<ul style="list-style-type: none">Remove the protection cap before use. Use this cap to protect shaft tubes from damage when not in service. Dispersing causes the medium to heat up. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. The function of the dispersion tools depends on the condition of the sharp edges of the rotor. Abrasive media can round off these edges fast, whereby the disperse effect diminishes.					
Designation	Application	Max. permissible speed [rpm]	Working range pressure [bar]	Working range vacuum [mbar]	
R 50	Wet crushing, suspension, emulsifying, deagglomerating (Batch-operation)	10,000	-	-	

Application instruction					
<ul style="list-style-type: none">Remove the protection cap before use. Use this cap to protect shaft tubes from damage when not in service. Dispersing causes the medium to heat up. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. The function of the dispersion tools depends on the condition of the sharp edges of the rotor. Abrasive media can round off these edges fast, whereby the disperse effect diminishes.					
Designation	Application	Max. permissible speed [rpm]	Working range pressure [bar]	Working range vacuum [mbar]	
R 50	Wet crushing, suspension, emulsifying, deagglomerating (Batch-operation)	10,000	-	-	

Designation	Application	Max. permissible speed [rpm]	Working range pressure [bar]	Working range vacuum [mbar]	
R 50	Wet crushing, suspension, emulsifying, deagglomerating (Batch-operation)	10,000	-	-	

Spare parts list					
Item	Designation	Item	Designation		
1	Protection cap	10	Shim		
3	Grooved pin	11	Rotary shaft		
4	Deep-groove ball bearing	13	Dissolver disc		
5	Circlip	14	Cage		
6	Circlip	16	Static ring		
7	Shim	17	Friction ring		
8	O-ring	18	O-ring		
9	Bearing bushing				

Spare parts diagram see front page.

For orders of spare parts please specify the fabrication number on the shaft tube, the type of dispersion tool and the designation of the spare part.

Item	Designation	Item	Designation		
1	Protection cap	10	Shim		
3	Grooved pin	11	Rotary shaft		
4	Deep-groove ball bearing	13	Dissolver disc		
5	Circlip	14	Cage		
6	Circlip	16	Static ring		
7	Shim	17	Friction ring		
8	O-ring	18	O-ring		
9	Bearing bushing				

Cleaning, disinfection and sterilization					
<p>For pre-cleaning the R 50 agitator shaft, operate it in in a detergent for dissolving the residuary of the substance. Due to the large flow rate, the rotor is cleaned to a large extent.</p> <p>The dissolver disc and cage can be removed for cleaning purposes.</p> <p>Hold the two flats with the open end wrench and unscrew the rotors with the other open end wrench. The protection cage can be removed by turning the screw with the allen wrench.</p> <p>Reassemble the dispersion tool in the reverse order.</p>					
Please note, that the protection cage is mounted correctly. The lower ring of the protection cage and the rotor must have the same level, see fig. 1.					

Technical Data						
Designation	Peripheral speed at 10,000 rpm	Min./max. immersion depth	Materials in contact with medium	Suitable for solvents	Temperature range upto max.	Vacuum/pressure upto
R 50	20,9-23,6 m/s	50-180 mm	AlSi 316 L	yes	100 °C	-

Warranty						
<p>In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.</p> <p>The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.</p>						

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