IKA

designed for scientists

IKA AOD 1

Betriebsanleitung	DE	3
Ursprungssprache		
Operating instructions	EN	16
Mode d'emploi	FR	29
使用说明	ZH	42

Indicaciones de seguridad	ES	55	Bezpečnostní upozornění	CS	69
Veiligheidsinstructies	NL	56	Biztonsági utasítások	HU	71
Norme di sicurezza	IT	58	Varnostna navodila	SL	72
Säkerhetsanvisningar	SV	59	Bezpečnostné pokyny	SK	73
Sikkerhedshenvisninger	DA	61	Ohutusjuhised	ET	75
Sikkerhetsinformasjon	NO	62	Drošības norādes	LV	76
Turvallisuusohjeet	FI	63	Saugos nurodymai	LT	77
Instruções de segurança	PT	65	Инструкции за безопасност	BG	79
Υποδείξεις ασφαλείας	EL	66	Indicații de siguranță	RO	80
Wskazówki bezpieczeństwa	PL	68	Указания по технике безопасности	RU	81

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Contents

	Page
1 User notes	17
1.1 EU Declaration of conformity	17
1.2 Explication of warning symbols	17
1.3 Notes on using these operating instructions	17
1.4 System features	17
1.5 Warranty	18
2 Safety instructions	18
3 Transport, Storage, Installation Location	20
3.1 Transport and storage conditions	20
3.2 Unpacking	20
3.3 Scope of supply	20
3.4 Installation location	20
4 Description of system components	21
4.1 Remote ignition unit AOD 1.2	21
4.2 Guard AOD 1.3	21
4.3 Ignition head	21
4.4 Further components	21
5 Setting-up and initial start-up	22
6 Carrying out decomposition tests	23
6.1 Recommendations for the sample	23
6.2 Use of combustible crucible C 14	23
6.3 Weighing the sample	23
6.4 Preparation of decomposition vessel AOD 1.1	24
6.5 Contact to and positioning the decomposition vessel	25
6.6 Decomposition	25
6.7 Sample transfer and absorption of combustion gases	25
6.8 Cleaning the decomposition vessel	26
7 Care and maintenance	26
7.1 General cleaning recommendations	26
8 Accessories and consumables	27
8.1 Accessories	27
8.2 Consumables	27
9 Technical data	27
9.1 Technical data for remote ignition unit AOD 1.2	27
9.2 Technical data for decomposition vessel AOD 1.1	27
9.3 Technical data for guard AOD 1.3	28
9.4 Technical data for oxygen filling station C 48	28
10 Index	28

1.1 EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/68/EU and 2011/65/EU and conforms with the following standards or normative documents: EN ISO 12100.

A copy of the complete EU Declaration of Conformity can be requested at sales@ika.com.

1.2 Explication of warning symbols

General hazard

This symbol identifies information **that is of absolute importance to ensure your health and safety**. Failure to observe this information may be detrimental to your health or may result in injuries.

This symbol identifies information **that is of important to ensure problem-free technical operation of the system**. Failure to observe this information may result in damage to the Decomposition System AOD 1.

This symbol identifies information that is important to ensure problem-free operation of decomposition tests and for working with the decomposition system. **Failure to observe this information may result in inaccurate measurement results.**

1.3 Notes on using these Operating Instructions

In this section you will learn how to make the most effective use of these Operating Instructions so as to be able to work safely with the Decomposition System AOD 1.

Section 7 contains important information on care and maintenance, and accessories, consumables, and technical data are given in Sections 8 and 9. The index is in section 10.

The instructions given in section 2 "Safety instructions" must be followed without fail!

Studying section 1-8

Work through sections 1 to 8 in numerical order. Section 3 "Transport, Storage, Installation Location" is particularly relevant to system reliability and ensuring high accuracy of measurements. Section 4 describes the components of the Decomposition System AOD 1.

Carrying out tests

Once you have carried out the procedures described in section 5 "Setting-up and initial start-up", the decomposition system AOD 1 is ready for trial decompositions. Section 6 "Carrying out decomposition trials" describes the complete test procedure from preparation of the components to the subsequent analysis.

In subsequent sections, the figures ①, ②, ③ etc. indicate instructions for operations which must always be carried out in the order given.

1.4. System features

Decomposition system AOD 1 is designed for the oxidative decomposition, underpressure, of solid and liquid samples that contain halogens and sulphur. It provides a fast, safe, low-maintenance method of quantitatively determining halogens and sulphur by decomposition.

The Decomposition system AOD 1 is built up from several components, whose function can also be extended with the aid of suitable accessories. To ensure the greatest possible safety during a decomposition test, **IKA** recommends use of the Guard AOD 1.3. You will find a description of the individual components in section 4 "Description of system components" or in their operating instructions.

The system has the following important features:

- In an atmosphere of pure oxygen, core temperatures of over 1300°C, at pressures up to 195 bar, can be reached in the crucible.
- The decomposition vessel is made from a highly halogen-resistant special alloy. For quantitative determination of fluorine, chlorine, bromine, iodine and sulphur, the inner surfaces and parts are catalytically activated.
- The gases generated during combustion can be dissolved in a special absorption solution, which is added to the decomposition vessel.
- Manual ignition of the sample using the remote ignition unit.
- Guard AOD 1.3 for safe operation (optional).

1.5 Warranty

In accordance with IKA Terms and Conditions of Sale, the warranty period is 12 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

- The detection method for dissolved ions can be selected by the user. Ion chromatography, detection with ion-selective electrodes, or titration are suggested methods. The following publications use these methods:
 - GIT 4/96: Determination of element specific halide- and sulfur-contents inorganic matrices.
 - GIT 7/96: New aspects within calorimetry.
 - GIT 11/96: Determination of calorific values with simultaneous halogen and sulfur decomposition.

These publications are available from **IKA**.

The Remote Ignition Unit AOD 1.2 must only be opened by an authorised service workshop. If servicing is required, we recommend that you take advantage of our customer service facilities. Further, please make yourself familiar with the relevant safety and accident prevention regulations.

Parts carrying electric voltage:

IKA accepts no liability for damage or costs that arise due to accidents, misuse of the unit, or unauthorised changes, repairs or modifications.

2

Safety instructions

Read the operating instructions in full before starting up and follow the safety instructions.

Keep the operating instructions in a place where they can be accessed by everyone.

Ensure that only trained staff work with the appliance.

Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Intended purpose:

The decomposition system AOD 1 is designed for the test sample decomposition of halogenated and sulphurous substances. The system must only be used for this purpose in conjunction with the original **IKA** reaction vessel AOD 1.1. For detailed information, see the operating manual for the reaction vessel.

Operating requirements:

The operator of this equipment must ensure safe operation of the decomposition vessel AOD 1.1 by installing a suitable safety device (e.g. safety guard AOD 1.3).

The maximum **energy input** in the decomposition vessel must not exceed **30000 J** and the sample mass should therefore be chosen accordingly. The maximum permitted **operating pressure** is **195 bar (19,5 MPa)**; do not exceed this pressure. The maximum permitted **operating temperature** is **50 °C**; do not exceed this temperature.

Do not overfill the decomposition vessel with the sample substance. Fill the decomposition vessel with oxygen at a pressure not exceeding the permitted maximum of **40 bar (4 MPa)**.

Check the pressure on the pressure reducer. Before the start of every combustion cycle, perform a seal test (following the instructions in the decomposition vessel operating manual).

When using protective device AOD 1.3, a **minimum distance of 2 metres** should be observed. Should a decomposition vessel burst, the device will not protect against damage to hearing. Always wear hearing protection in order to avoid damage to the ears.

Explosivstoffe:

Some substances tend to combust in an explosive manner due to peroxide build-up. This can cause the decomposition vessel to burst.

The decomposition vessel AOD 1.1 must not be used for the testing of explosive samples.

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Substances whose combustion behaviour is unknown must be tested prior to combustion in the decomposition vessel AOD 1.1 to establish their combustion behaviour (**explosion hazard**). When you burn an unknown sample, maintain a safe distance from the decomposition vessel. With the guard AOD 1.3 fitted, the minimum safe distance is 2 metres. Benzoic acid must only be burnt in its compressed form. Combustible dusts and powders must be pressed before testing. Explosion hazard: over-dried dusts and powders (e.g. wood shavings and sawdust, hay and straw) burn in an explosive manner. Wet these substances before testing. Keep slightly flammable fluids with low vapour pressure (tetramethyl-dihydrogendisiloxan) away from cotton yarn; prevent all contact between these two substances.

Observe the relevant health and safety regulations. Always wear personal protection equipment (PPE).

Combustion residue, auxiliary materials:

The inside wall of the decomposition vessel may become coated with toxic combustion residues such as gas, ash and other deposits.

When handling combustion samples, combustion residues and auxiliary materials, the appropriate safety requirements must be observed. The following are examples of substances that may cause dangers: corrosive, easily flammable, capable of exploding, contaminated with bacteria, toxic.

Oxygen:

When working with oxygen, observe the appropriate requirements. **Danger warning:** As a compressed gas, oxygen promotes combustion, supports combustion intensively and may react violently with combustible substances. **Do not use any oil or grease!**

Using a crucible made of stainless steel:

When using crucibles made of stainless steel, their condition should be carefully checked after every experiment. A reduction in the thickness of the material may cause the crucible to burn and may damage the decomposition vessel.

For reasons of safety, crucibles must not be used any more after a maximum of 25 combustion procedures.

Ignition wire:

The decomposition vessels are equipped with a fixed ignition wire. Since the ignition wire is subjected to a high degree of material stress during combustion, it should be replaced after 50 tests at the latest.

Specification of the decomposition vessel:

The decomposition vessel has been manufactured in accordance with the European Pressure Equipment Directive 97/23/EC. The decomposition vessel has been **pressure tested to 280 bar (28 MPa)** and seal tested with oxygen to 30 bar (3 MPa).

The decomposition vessel is a laboratory **autoclave** and must be tested by a skilled person after each single use. Single use here is understood also as a series of tests where approximately the same pressures and temperatures are used. Laboratory autoclaves must only be operated in special chambers or behind protective walls.

Regular test intervals:

The decomposition vessel must be subject to inner tests and pressure tests performed by a **skilled person** at regular intervals. The test intervals are to be defined by the user on the basis of experience, operating conditions and the type of samples tested.

The declaration of conformity on this equipment becomes void if mechanical modifications are made to the laboratory autoclave or where the resistance of the vessel can no longer be guaranteed because of very severe corrosion. An example of very severe corrosion is pitting caused by halogen.

The **thread** on the body of the decomposition vessel and the union nut are subject to high stresses and must therefore be checked for **wear** at regular intervals.

The condition and the efficiency of the seal must be tested using a **seal test** (see the operating manual for the decomposition vessel). Pressure testing and servicing must be performed by **skilled personnel** only.

The manufacturer hereby specifies that the decomposition vessel must be tested or returned to our works for repair, after 1,000 test hours, after one year or earlier if the operating conditions warrant this.

Definition of person with professional training:

A person with professional training as defined in these operating instructions is someone:

1. whose training, knowledge and experience gained through practical activities ensures that that person will perform the tests in a proper manner.
2. who is sufficiently reliable.
3. who is not subject to any instructions in terms of testing activity.
4. who is equipped with suitable testing equipment if necessary.
5. who can provide suitable proof demonstrating compliance with the requirements listed in 1.

Operating pressure containers:

The operation of pressure containers is regulated by national directives, standards and regulations. Operators of pressure vessels are required to keep these vessels in good condition, operate them correctly, test them correctly, carry out the necessary servicing and maintenance and to take the necessary safety precautions.

A pressure vessel must not be operated when it is faulty or where it can cause hazard to the operator or third parties.

The European Pressure Equipment Directive is available from the publisher Beuth Verlag.

Use only original spare parts from IKA!

3 Transport, Storage, Installation Location

3.1 Transport and Storage Conditions

During transport and storage, the system must be protected against mechanical shock, vibration, dust deposits, and corrosive atmospheres. In addition, the relative humidity should not exceed 80%. Transport should only be undertaken with the unit in its original packing.

3.2 Unpacking

Please unpack system components carefully and check for any signs of damage. It is important that any transport damage is noted during unpacking. If necessary, the damage must be assessed immediately by the transport company (post, railway or transport contractor).


3.3 Scope of Supply

The standard scope of supply of the decomposition system AOD 1 consists of:

- 1 x Remote ignition unit AOD 1.2 (see item 1 in illustration) with ventilation grip,
- 1 x Decomposition vessel AOD 1.1 (see item 2 in illustration) with fixed platinum ignition wire C 5012.3 and a carrying grip,
- 1 x Oxygen filling station C 48 (see item 4 in illustration),
- 1 x IKA control standard for chlorine and sulphur AOD 1.11 (not illustrated),
- Operating instructions for decomposition system AOD 1, oxygen filling station C 48 and decomposition vessel AOD 1.1 (not illustrated).

The decomposition systems AOD 1 can be extended as required with the following optional accessories:

- Guard AOD 1.3 with ignition cable connected, cable length 5 m (recommended, see item 3 in illustration),
- Remote ignition head AOD 1.13 with ignition cable connected, cable length 5 m (is required when Guard AOD 1.3 is not used),
- Venting station C 7030 with gas-washing bottle to DIN 12596 for gas absorption (see item 5 in illustration),
- Reducing valve C 29 (not illustrated),
- Pelleting press C 21.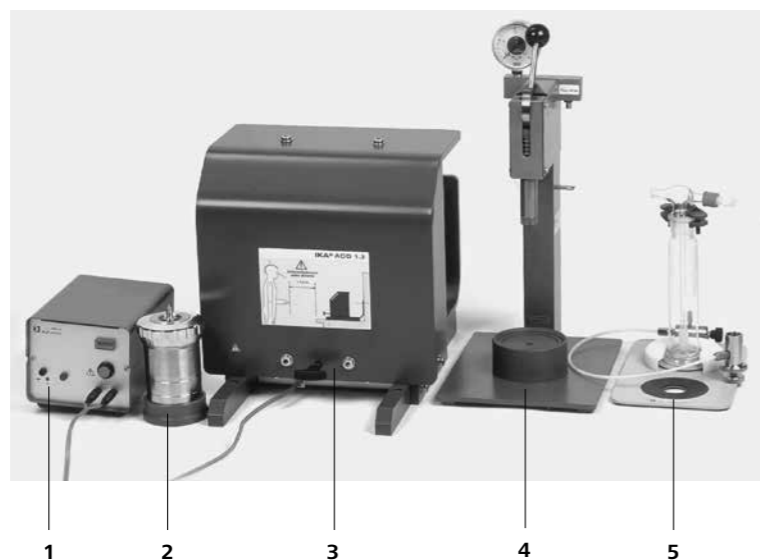

3.4 Installation Location

When installing the AOD 1 System, observe the national and local regulations for the operation of pressure vessels that apply at the site selected. The site selected for installation of the AOD 1 System must be chosen so that the guard screens people from the decomposition vessel during a test. The responsibility for safe operation of the system always rests with the operating authority. Please observe Section 1 "Safety instructions".


Recommendation for operation with Guard AOD 1.3:
The Guard AOD 1.3 must be positioned so that no one can stay in the unprotected zone behind it.

Recommendation for operation with another type of guard:
The operator must ensure that the site selected for installation of the system permits safe operation, particularly taking into account features of the guard being used.

The AOD 1 system must be set up on a bench or in a fume cupboard to ensure clean working conditions. The system requires an electrical supply corresponding to the data on the type plate of the remote ignition unit and an oxygen supply (99.95 % pure oxygen, quality 3.5; pressure 30 bar) with a pressure gauge where it is installed. The oxygen supply must comply with local regulations. Please read the operating instructions for the oxygen filling station C 48. There must be a shut-off valve for the oxygen supply. Follow the instructions for oxygen in Section 1 "Safety instructions".



4 Description of System Components

4.1 Remote ignition unit AOD 1.2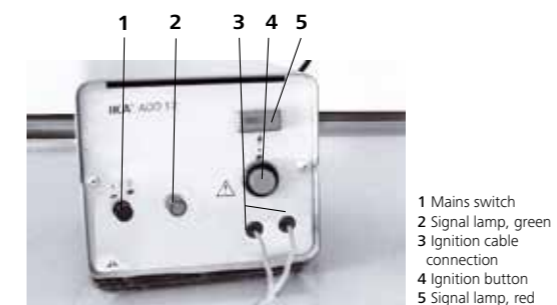

The remote ignition unit provides the current necessary to ignite a test sample. After the ignition button has been pressed, the ignition wire in the decomposition vessel is heated so that the cotton thread or the combustible crucible, and thus the sample, are ignited.

4.2 Guard AOD 1.3

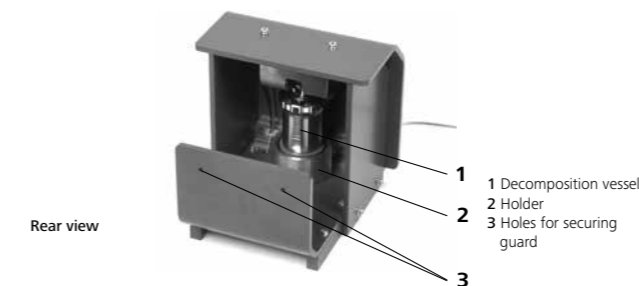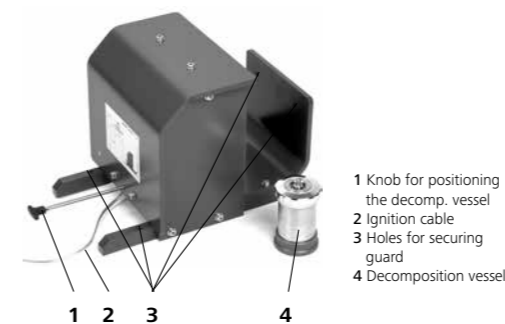The Guard AOD 1.3 with its integrated ignition contact helps to protect persons working with decomposition vessels. However, protection is only provided for persons who are in front of the guard. When the Guard AOD 1.3 is used, a sample can only be ignited when the decomposition vessel is correctly positioned within the guard. The decomposition vessel must be placed in the holder, and pulled to the ignition position using the knob.

4.3 Ignition head

The Ignition head is required if Guard AOD 1.3 is not employed. It has an ignition cable that is connected to the remote ignition unit. Placing the ignition head on the decomposition vessel makes an electrical connection to the ignition wire in the vessel.

4.4 Further Components

Information on the decomposition vessel AOD 1.1, the oxygen filling station C 48 and the venting station C 7030 are given in their operating instructions.

First the components of the decomposition systems AOD 1 must be unpacked and brought to the selected location. This location must comply with the requirements for safe operation described in section 3.4 "Installation Location". In addition, distilled water, dilute nitric acid for cleaning purposes, 0.25-molar caustic soda, and 30% hydrogen peroxide, all in high-purity form, should be readily available. Carry out the following steps:

①

Connection of the Oxygen Filling Station C 48

The oxygen pressure should be at least 30 bar, but must not exceed 40 bar. Quality 3.5 oxygen (99.95 % pure) must be used.


Connect the Oxygen Filling Station to your laboratory oxygen supply using the hose provided. Further details are given in the operating instructions for the Oxygen Filling Station C 48.

②

Positioning the guard

• Operation with Guard AOD 1.3

Place the Guard AOD 1.3 with its opening to the rear at the intended location, and secure it there so that it cannot be displaced. To this end the Guard has a number of holes in its supports and rear surface for screws or other fasteners.

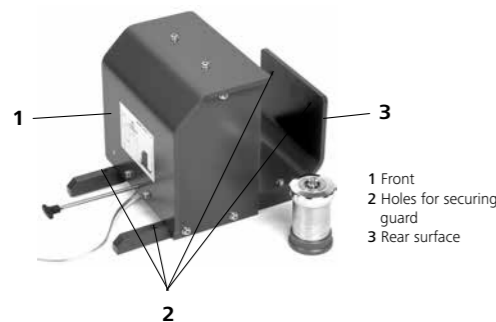• Operation with another guard

Place the guard at the intended location, and secure it there so that it cannot be displaced. The guard must be strong enough to prevent injury to persons if the decomposition vessel bursts.

③

Connecting the ignition cable to the remote ignition unit

The plugs on the ignition cable must only be connected to Remote Ignition Unit AOD 1.2.

• Operation with Guard AOD 1.3

The guard AOD 1.3 has an ignition cable 5 m long. Connect guard AOD 1.3 with the remote ignition unit AOD 1.2 by connecting the two plugs on the cable to the sockets provided on the remote ignition unit.

• Operation with another guard

Connect the ignition head to the remote ignition unit AOD 1.2 by connecting the two plugs on the ignition cable to the sockets provided on the remote ignition unit.

④

Setting-up, connecting, and switching on the Remote Ignition Unit

Position the remote ignition unit so that while you are operating it you are protected by the guard. The remote ignition unit must be at least 2 metres from the guard. Check that the voltage on the type plate of the remote ignition unit corresponds to the mains supply you will be using. If the supply is correct, connect the mains lead to it. Switch the remote ignition unit on at the mains switch. When the unit is switched on, the green signal lamp illuminates.

⑤

Initial use of the decomposition vessel AOD 1.1

The decomposition vessel is supplied closed. Before using it for the first time, un-screw the ring nut and use the grip to remove the lid of the decomposition vessel (see also the operating instructions for decomposition vessel AOD 1.1). Clean all parts of the decomposition vessel with dilute nitric acid, and rinse them in distilled water until all traces of acid have been removed.

Internal parts and the internal surfaces have a cloudy and partly stained appearance. This condition is of great importance for the subsequent analysis. It indicates catalytically-active surfaces, and must not be removed by brushing or using hard tools. Wipe out and dry the decomposition vessel only with a soft, lint-free cloth. Further information is given in the operating instructions for decomposition vessel AOD 1.1. Clean the active parts of the decomposition vessel before first using it by carrying out two combustion runs, each with two benzoic acid tablets (1 g), as described in the next section "Carrying out decomposition trials".

The prerequisites for reproducible and accurate testing are cleanliness when handling the decomposition vessel, purity of the chemicals used, and meticulously following the chosen procedure. Always observe the hints and instructions in section 2 "Safety instructions" and the following sections; they are there to ensure safe operation.

6.1 Recommendations for the sample

Samples to be analysed must be homogenised or, if necessary, ground before they are weighed. Samples that readily generate dust are easier to handle if they are first pressed into pellets. This also helps achieve more uniform combustion.

Please also read the safety notes for decomposition vessel AOD 1.1.

Solids

Normally, solid fuels can be burnt directly in powder form. Fast-burning substances (e.g. benzoic acid) must not be burnt in a loose form.

Fast-burning substances tend to spluttering combustion. Complete combustion is then no longer certain. In addition, the inner wall of the decomposition vessel may be damaged. Such substances must be pressed to form pellets before they are burnt.

The IKA C 21 Pelleting Press is suitable for this task (accessory).

Liquids

Most liquids can be weighed directly into the crucible. Liquids that are cloudy or with separable water must be dried or homogenised before weighing. For solvents with different phases, each phase should be tested separately.

Highly-volatile substances

Loss due to evaporation, dusting, and environmental factors can be minimised by weighing the sample in acetobutyrate C 10 or gelatine capsule C 9 (see Accessories). These capsules also act as a combustion aid.

For a low-flammability substance, a combustion aid should be weighed into the crucible together with the sample.


Combustion aids

Low-calorific-value samples may also require an aid to combustion such as paraffin or benzoic acid for thorough burning. Paraffin can also be used to treat samples that tend to produce dust, acting as a binding agent and improving combustion. Besides the capsules mentioned above, polyethylene combustion bags C 12 and C 12A or combustible crucible C 14 (accessories) can also be used.

At low detection levels, it is better to avoid using a cotton thread as an ignition aid, and to use a paraffin strip C 15 instead. In this way, the effects of impurities in the cotton, which adversely affect the lower limits of detection, can be minimised. We strongly recommend carrying out blind determinations to establish whether impurities are likely to be present.

6.2 Use of Combustible Crucible C 14

The decomposition vessel AOD 1.1 provides for the use of combustible crucibles. Using combustible crucible C 14 considerably improves the burning of a sample. The combustible crucible is made of acetobutyrate and has an energy input of 5000 Joule in the decomposition vessel. When a combustible crucible is used, a cotton thread is not required.

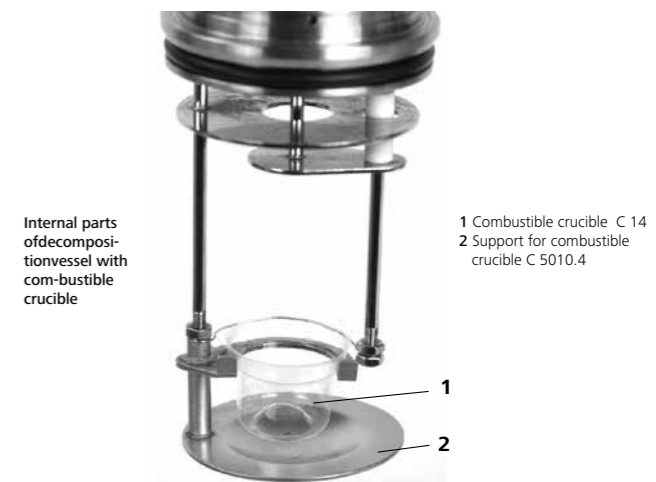When a combustible crucible C 14 is used, the decomposition vessel must have the Support for combustible crucible C 5010.4 (accessory). The sample can then be weighed directly into the crucible. The combustible crucible must be placed in the crucible holder so that it is in direct contact with the ignition wire. To ensure this, clip the combustible crucible under the ignition wire.

When ignition is initiated and the wire starts to glow, the crucible is ignited, and the sample burnt from all sides. The crucible is completely destroyed. Experience has shown, that the additional energy input from a combustible crucible, with the supply of oxygen from all sides, results in better combustion than in a quartz crucible.

The combustible crucible has the same dimensions as a quartz crucible C 4. Even if the AOD decomposition vessel is equipped for combustible crucibles, a quartz crucible can still be used.

6.3 Weighing the Sample

Before every decomposition trial, the crucible and decomposition vessel must be thoroughly rinsed in distilled water and then dried. When weighing the sample use only clean implements. Avoid all skin contact with internal parts of the system – this can be a cause of false results.

① **Sample weighing**

The operating temperature of the decomposition vessel must not exceed 50 °C. This corresponds to a maximum energy input of ca. 20000 J. Choose the sample mass accordingly. If the energy input is too large, the decomposition vessel may be damaged.

With damaged decomposition vessels, there is a danger of bursting! Always observe the operating instructions for the decomposition vessel!

When working with an unknown substance, start with a very small sample to determine its energy potential, and keep a safe distance from the guard.

The weight of the sample will depend on the test requirements and the halogen or sulphur content of the substance – it may vary between a few milligrams and up to 1 g.

6.4 Preparation of Decomposition Vessel AOD 1.1

 Decomposition Vessel AOD 1.1 is supplied with a fixed platinum ignition wire C 5012.3. This fixed ignition wire is, however, subject to wear, and should be checked after every combustion test. If the wire has become visibly thinner, it must be replaced (see the operating instructions for Decomposition Vessel AOD 1.1).

① **Solution in decomposition vessel**

Ten ml of 0.25 molar NaOH solution and 100 to 200 µl of 30 % hydrogen peroxide solution are standard additions to the decomposition vessel. Other additions and concentrations are also possible, depending on the anticipated composition of the sample (see also instructions for using AOD 1.11 and AOD 1.12).

② **Securing a sample in the holder**

Then a crucible with the weighed sample is placed in the crucible holder, and secured with the retaining ring. Place the retaining ring below the crucible holder around the crucible. The crucible is then held to the holder, and cannot come loose, even when the vessel is shaken.

 If a combustible crucible is used, the retaining ring is not required.

③ **Closing the decomposition vessel**

Close the decomposition vessel by placing the lid with the sample in it until there is metallic contact. Screw the ring nut on to the decomposition vessel until it is just tight. The decomposition vessel is now ready for filling with oxygen.


④ **Filling the decomposition vessel using Oxygen Filling Station C 48**

Fill the decomposition vessel with oxygen using the Oxygen Filling Station C 48. Please read the C 48 operating instructions.

The oxygen pressure should be at least 30 bar, but must not exceed 40 bar. Quality 3.5 oxygen (99.95 % pure) should be used.

Hold the hand lever of the Oxygen Filling Station firmly while moving it; the pressure in the decomposition vessel can cause the lever to move rapidly upwards (danger of shock).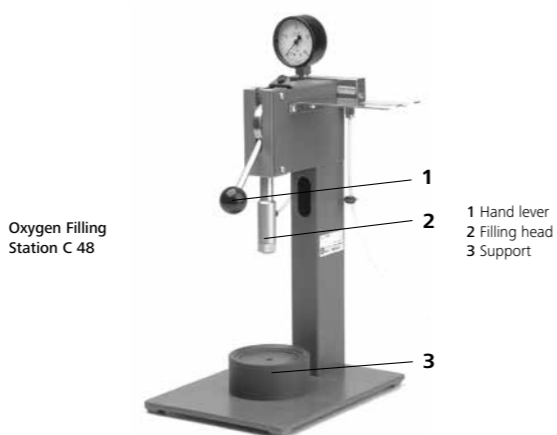

Place the decomposition vessel that is to be filled on the support. Operate the hand lever to move the filling head down to the decomposition vessel, and push it over the filling nozzle. Take care that the decomposition vessel is correctly positioned. If the vessel is not correctly positioned, there will be a loud noise of escaping gas during the filling operation. There will be a similar noise if the decomposition vessel leaks due to wear of the seals (see Section 7 "Care and Maintenance").

After about 40 seconds, the decomposition vessel will have been filled, and can be removed from the oxygen filling station. To do this, move the hand lever upwards; the oxygen supply is cut off and filling stops.

The decomposition vessel is now prepared for carrying out a test.

6.5 Contact to and positioning the decomposition vessel

Make sure that the guard has been installed as described in Section 3.4 "Installation Location" and Section 5 "Setting-up and Initial Start-up".

Operation with guard AOD 1.3

Place the decomposition vessel behind the Guard AOD 1.3 in the holder (see Section 4.2 "Guard AOD 1.3" for illustration). With the knob on the front of the guard, draw the decomposition vessel into the guard until the stop is reached at the ignition position. In this position, the decomposition vessel is in contact with the ignition contact of the guard and thus – as long as the ignition cable is connected to the Remote Ignition Unit as described in Section 6 – with the remote ignition unit.

Operation with another guard

When positioning the decomposition vessel behind the guard, make sure that no one can be injured by flying debris if a decomposition vessel bursts.

The ignition head (see Section 4.3 "Ignition Head") is placed on the decomposition vessel and locked by turning it anti-clockwise. This makes the electrical connection to the remote ignition unit.

Then position the decomposition vessel behind the guard you are going to use.

6.6 Decomposition

Before initiating decomposition at the remote ignition unit, during decomposition, and right through to the end of the cooling phase, make sure that all persons (including all their limbs) are within the area protected by the guard. Maintain a safety distance of 2 m to the guard.

For removal and transfer of the decomposition vessel to a water bath, use the carrying grip included in the scope of supply. Direct contact with the decomposition vessel after a test can cause burns.

To ignite a sample, press the red button (see illustration in Section 4.1 "Remote Ignition Unit") on the remote ignition unit once. A warning signal sounds.

Following remote ignition, wait at least one minute before removing the decomposition vessel again. The combustion process will heat up the decomposition vessel. At the end of the waiting period, lock the carrying grip on to the lid of the decomposition vessel by turning it anti-clockwise. You can now use the carrying grip to transfer the decomposition vessel to a water bath. For cooling and quantitative transfer of the reaction gases into solution, the decomposition vessel should be completely submerged in the water bath for about 5 minutes.

 If the red signal lamp illuminates during ignition, the ignition wire is defective. If the current is too high, the fuse on the rear of the unit will trip out. After about one minute, the fuse can be reset by gentle finger pressure. The unit is then ready for use again.

6.7 Sample transfer and absorption of combustion gases

When venting the decomposition vessel, make sure that combustion gases cannot escape into the room.

On completion of a test, the decomposition vessel must be vented. Venting can be carried out using the ventilation grip, which is included in the scope of supply, or with the Venting Station C 7030 (accessory). Depending on the sample composition, venting the combustion gases without absorption can yield low values for halogen and sulphur content. For quantitative decomposition, it is therefore essential to pass the combustion gases through an absorption solution. In these cases, venting should be carried out using Venting Station C 7030.

① **Venting the decomposition vessel**

While the decomposition vessel is still under pressure, it should be shaken gently to aid absorption of the residual gases; a homogeneous distribution of the substance to be analysed and the condensate in the liquid phase is a prerequisite for the subsequent analysis.

• **Venting with the venting grip:**

Place the venting grip on the decomposition vessel and lock it by turning it anticlockwise. Vent the vessel in a fume cupboard by pressing the knob.

• **Venting with Venting Station C 7030:**

Venting using the Venting Station C 7030

To vent a decomposition vessel using the Venting Station C 7030, please follow the operating instructions for the Venting Station C 7030.

 Slow venting through a washing bottle is, for example, essential for iodine determination.

② **Checking for complete combustion**

Open the decomposition vessel and check the crucible and decomposition vessel wall for signs of incomplete combustion. If you find evidence of incomplete combustion, reject the results of that test, and repeat it.

 Incomplete combustion can be caused if the oxygen pressure in the decomposition vessel is too low, or it indicates that a combustion aid is required.

③

Preparation for analysis

Transfer the absorption solution carefully to a measuring flask with distilled water. All components of the decomposition vessel must be thoroughly rinsed. For analysis with ion chromatography, it is recommended that dissolved carbon dioxide is removed from the absorption solution.

Alternatively, a quick but less accurate determination can be carried out by filling a decomposition vessel with a measured quantity of distilled water to obtain a "defined volume" (e.g. 20 ml or 100 ml). This can then be subjected to gravimetric or volumetric analysis.

After the distilled water has been added, the vessel must be closed again, and, with the aid of the Oxygen Filling Station C 48, filled with oxygen to about 3 bar. This is necessary to close the valve on the vessel. The vessel must then be vigorously shaken. A homogeneous distribution of the substance to be analysed and the condensate in the liquid phase is a prerequisite for the subsequent analysis.

After venting again, the ions and elements of interest can be determined using a suitable method of analysis.

④

Analysis

The detection method for dissolved ions can be selected by the user. Ion chromatography, detection with ion-selective electrodes, or titration are suggested methods. Particularly when the samples being tested contain other halogens (iodine, fluorine or bromine) besides chlorine and sulphur, and very consistent detection is required, ion chromatography is to be preferred.

For further information, please refer, for example, to DIN 38414 Part 18, Determination of AOX in sludges and sediments". IKA can provide further information about Hg and As determination using AAS spectroscopy.

6.8 Cleaning the decomposition vessel

If there is a suspicion that a sample, or the residues or gaseous products of combustion could present a health hazard, then protective clothing (e.g. gloves, breathing mask) must be worn. Combustion residues which are a hazard to health or the environment must be disposed of as hazardous waste. We refer you explicitly to the valid regulations.

For accurate results, it is absolutely essential that the decomposition vessel is clean and dry. After every test, the inner walls of the vessel, the internal fittings (mountings, electrodes etc.) and the combustion crucible (internally and externally!) must be thoroughly cleaned.

Inner walls of vessel The inner walls of the vessel and internal parts should be cleaned with distilled water or dilute nitric acid, and then carefully wiped out with an absorbent lint-free cloth. If the decomposition vessel cannot be properly cleaned using the measures described (e.g. due to burning, pitting, corrosion etc.), do not attempt to clean it mechanically. In such cases, contact our Technical Service!

Crucible Combustion residues in the crucible, e.g. carbon or ash, must also be wiped off carefully with an absorbent, lint-free cloth.

7**Care and Maintenance**

If the decomposition system is handled properly, only the seals of the decomposition vessel, oxygen filling station, and venting station require maintenance. All these units have O-ring seals. These are subject to wear and must be replaced when leakage occurs. The necessary O-rings are available in sets for each component.

Before every decomposition test, it is absolutely essential to make a visual inspection of all parts of the decomposition vessel. If there are signs of corrosion, mechanical damage, loose electrodes, or consumption of the ignition wire, tests must not be carried out.

For maintenance of the decomposition vessel, see the AOD 1.1 operating instructions!

7.1 General cleaning recommendations

- For cleaning disconnect the mains plug!
- Use only cleaning agents which have been approved by IKA to clean the devices:
Water containing surfactant / isopropyl alcohol.
- Wear protective gloves during cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Notes:

Electrical equipment must not be immersed in a cleaning agent.

We recommend the wearing of protective gloves during cleaning.

The operating authority is responsible for ensuring that appropriate decontamination measures are taken if hazardous materials are spilt on or in the instrument.

Before using any method for cleaning or decontamination other than those recommended by the manufacturer, consult the manufacturer to make sure that the intended method will not damage the instrument.

When replacing the mains cable, use a product of equivalent quality and performance.

Spare parts order

When ordering spare parts, please give:

- device type,
- serial number, see type plate,
- position number and description of spare part, see www.ika.com.

Repair

Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.

For repair, please request the "Safety Declaration (Decontamination Certificate)" from IKA or use the download printout of it from IKA website at www.ika.com.

If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

8**Accessories and consumables****8.1 Accessories****Ordering description**

AOD 1.3	IKA Guard
AOD 1.13	Remote Ignition Head (required when Guard AOD 1.3 is not used)
C 21	Pelleting press
C 29	Pressure reducer
C 5010.4	Support for combustible crucible
C 7030	Venting station with gas washing bottle to DIN 12596 for gas absorption

8.2 Consumables**Ordering description**

AOD 1.11	IKA control standard for chlorine and sulphur
AOD 1.12	IKA control standard for fluorine and bromine
C 710.4	Cotton threads, cut to length (500 pieces)
C 5012.3	Platinum ignition wire, replacement (2 pieces)
C 4	Quartz dish
C 9	Gelatine capsules
C 10	Acetobutyrate capsules (100 pieces)
C 12	Combustion bags, 40 x 35 mm (100 pieces)
C 12A	Combustion bags, 70 x 40 mm (100 pieces)
C 43	Benzoic acid (NBS 39i, 30 g)
C 43A	Benzoic acid (100 g)
C 723	Benzoic acid tablets (50 pieces)
C 14	Combustible crucibles (100 pieces)
C 15	Paraffin strips (600 pieces)

9**Technical Data****9.1 Technical data for remote ignition unit AOD 1.2**

Rated voltage/frequency	V	115, 50/60 Hz
	V	230, 50/60 Hz
Power consumption	W	45
Fuse	A	0,3 (automatic)
Degree of protection to DIN 40 050		IP 21
Protection class		1 (with protective earth)
Over-voltage category		2
Contamination level		II
Ambient temperature	°C	5... 40
Max. ambient relative humidity	%	80
Dimensions (B x D x H)	mm	135 x 185 x 115
Weight	kg	2,7
Housing		Sheet metal, painted

9.2 Technical data for decomposition vessel AOD 1.1

Please see the Operating Instructions AOD 1.1 for technical data of the decomposition vessel.

9.3 Technical data for guard AOD 1.3

Dimensions (B x D x H)	mm	300 x 530 x 320
Wall thickness	mm	10
Weight	kg	12,2
Length of connection lead	m	5 (only connection to remote ignition unit AOD 1.2)
Application		suitable for all Type AOD 1.1 decomposition vessels

9.4 Technical data for oxygen filling station C 48

Please see the Operating Instructions for the Oxygen Filling Station C 48.

Subject to technical changes!

10

Index

A

Absorption solution.....2-2, 6-6
Acetobutyrate capsule..... 6-1
Analysis2-2, 6-7
Analysis methods 6-7

C

Cloudy liquids 6-1
Combustible crucible 6-2
Combustion aid 6-1
Combustion bags 6-1
Consumption of ignition wire 7-1
Corrosion..... 7-1
Cotton thread.....6-1, 6-2

D

Detection method2-2, 6-7

E

Explosive substances..... 1-1

F

Fast-burning substances 6-1
Features 2-2
Fuse 6-6

G

Gelatine capsule..... 6-1

I

Ignition position4-2, 6-5
Impurities..... 6-1
Incomplete combustion..... 6-7
Intended purpose 1-1
Iodine determination..... 6-7
Ion chromatography.....2-2, 6-7

L

Leakage.....6-5, 7-1
Liquids6-1
Loose electrodes7-1

M

Mechanical damage.....7-1

O

Oxygen supply3-2

P

Paraffin strip.....6-1
Pressure containers.....1-3
Pressure vessels3-2

S

Safe operation 1-1, 3-2, 6-1
Sample6-3
Solids.....6-1
Standard additions.....6-4
Switching on5-2

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited
IKA Works (Thailand) Co. Ltd.