IKA

designed for scientists

IKA C 25

BETRIEBSANLEITUNG	DE	3
MANUAL INSTRUCTIONS	EN	4
MODE D'EMPLOI	FR	5
INSTRUCCIONES DE MANEJO	ES	6

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Warranty

In accordance with IKA Terms and Conditions of Sale, the warranty period is 12 months. For claims under the warranty please contact your local dealer. You may also send the device direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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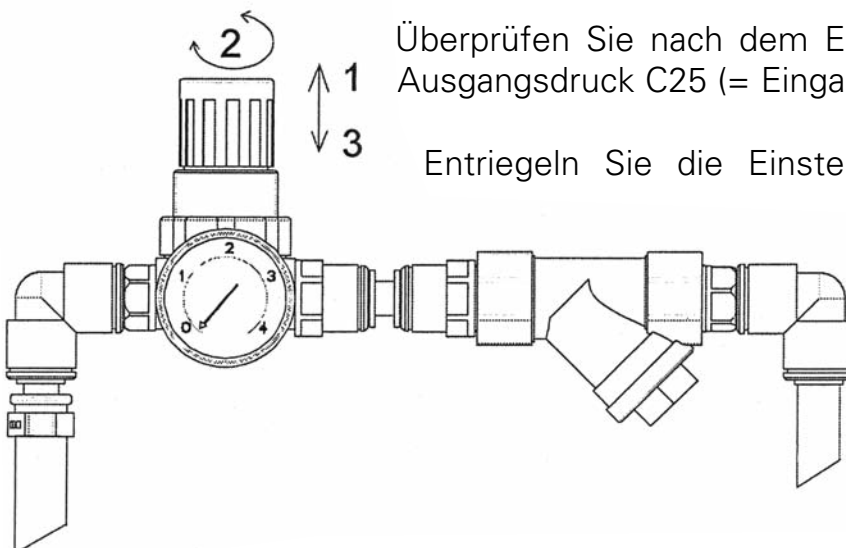
The IKA C25 Pressure Control Valve is a membrane pressure controller for laboratory devices that require a defined water input pressure (for example the IKA C2000 basic / control calorimeter).

C25 input pressure	max. 20 bar
C25 output pressure	0,5 - 3 bar
Operating medium	Water
Installation location	Any
Ambient temperature of medium	max. 50 °C
Connection	Ø 10 (G 1/4)

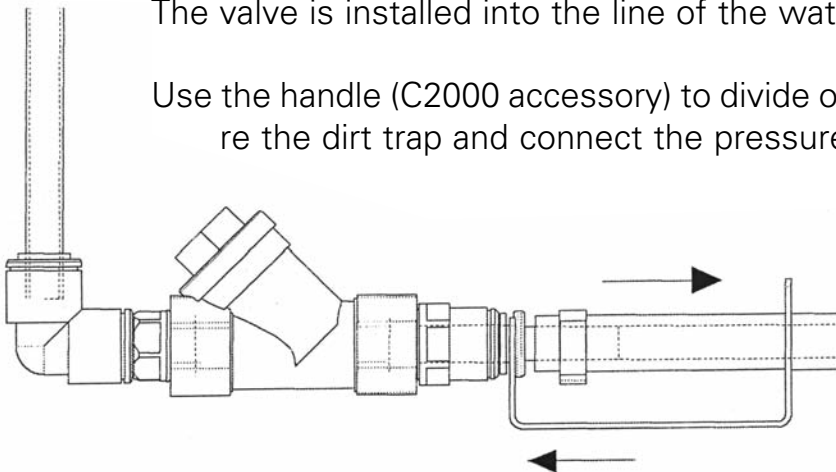
Commissioning

Commissioning with the IKA C2000 basic / control calorimeter

The C25 pressure valve is absolutely essential on the water faucet for operation of the C2000, and it must be preset to an output pressure of approximately 1,5 bar.


The valve is installed into the line of the water connection.

Use the handle (C2000 accessory) to divide off the hose of the water line before the dirt trap and connect the pressure control valve with the dirt trap.

Connect the hose onto the pressure control valve.

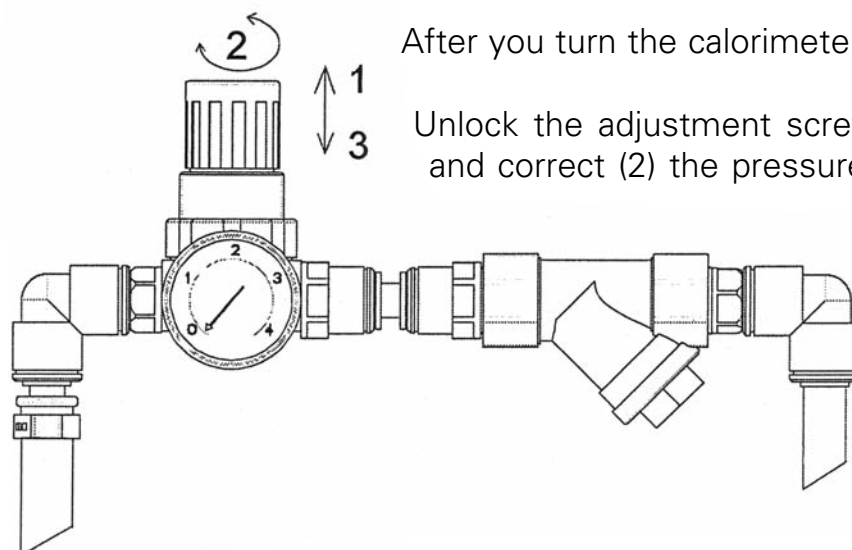The connections are set up correctly if you overcome a distinct resistance and plug the line in all the way to the stopper.

After you turn the calorimeter on, check the pre-pressure.

Unlock the adjustment screw by raising the hand wheel (1) and correct (2) the pressure to 1 - 1,5 bar if necessary (see the operating instructions for the C2000 basic /control).

Lock the hand wheel (3) in place so that it cannot be adjusted unintentionally.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited
IKA Works (Thailand) Co. Ltd.