D

**Installationsanleitung
Akku**

Seite 2

GB

**Installation Information
Rechargeable battery**

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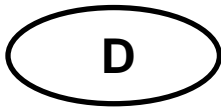
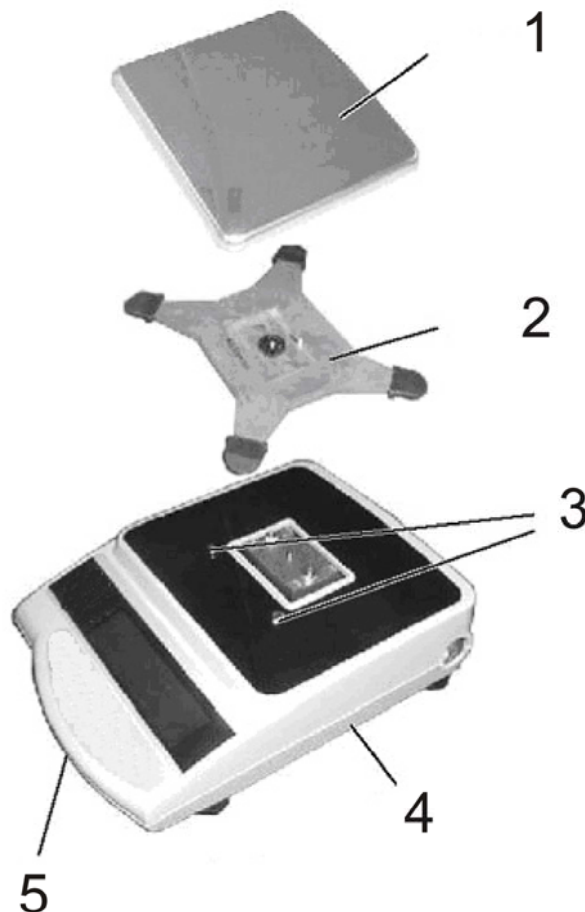
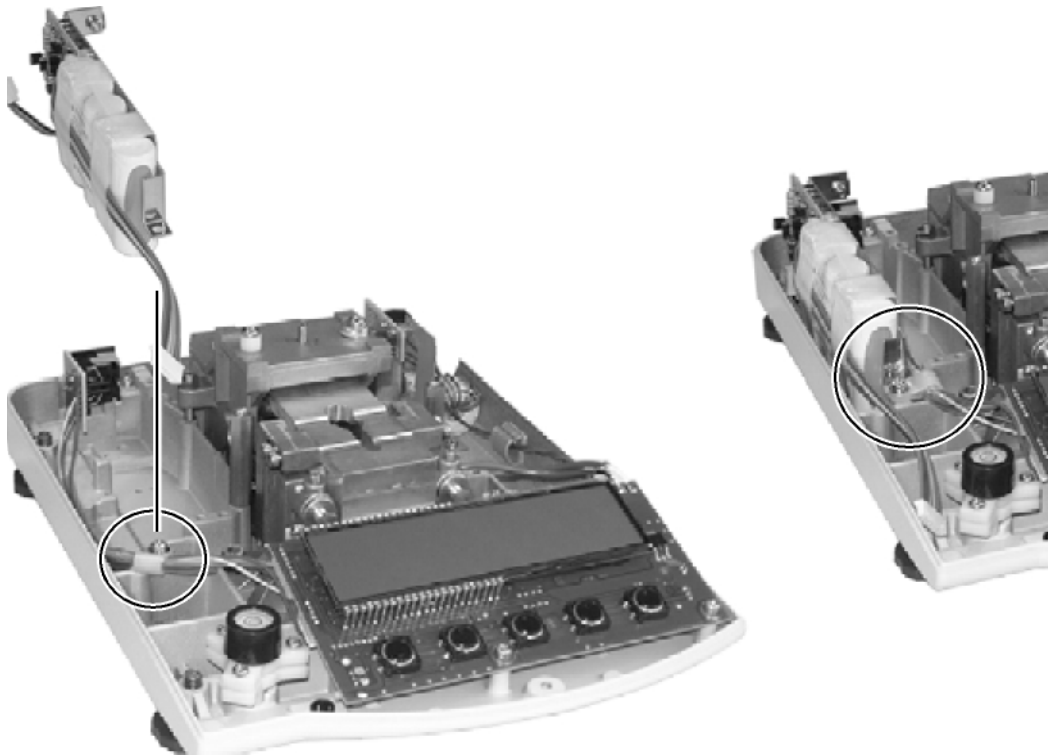
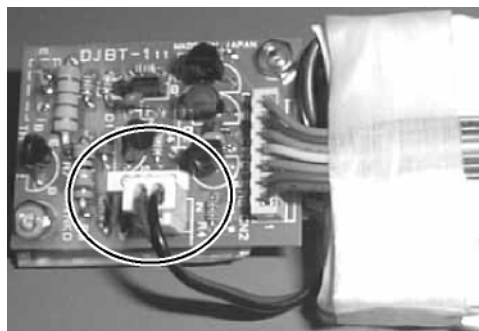
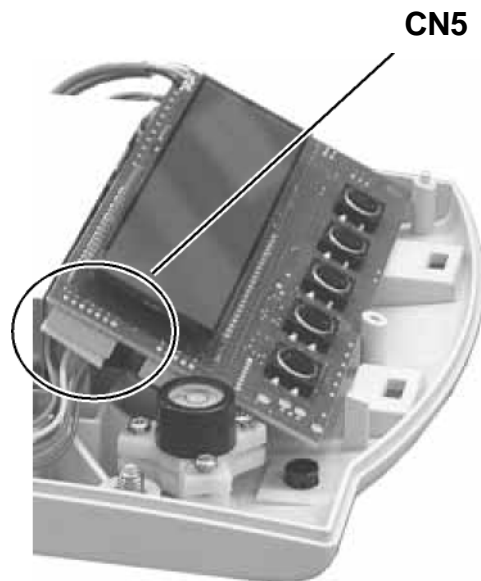
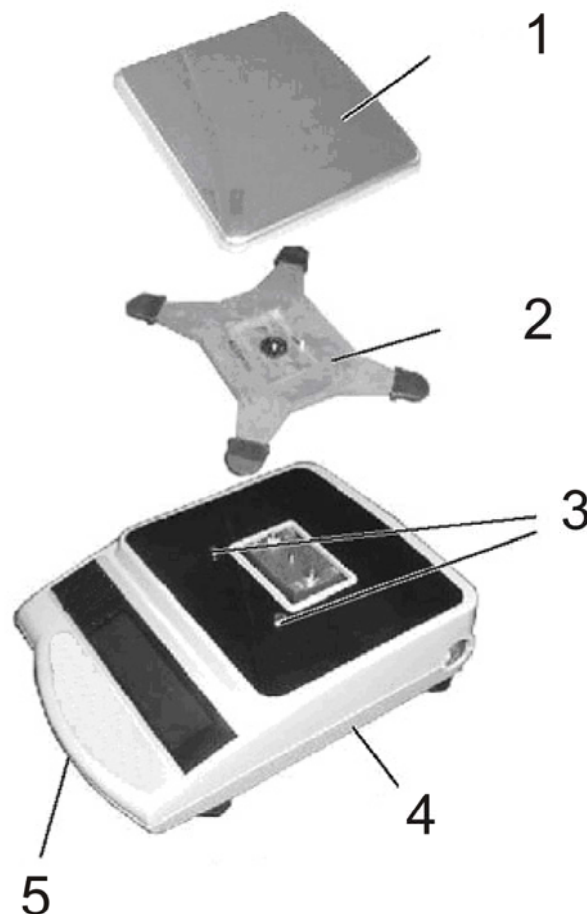
KERN EG-A06 for models with max ≥ 2200 g

How to remove the case:


- Disconnect the balance from the mains
- Remove weighing plate (1) and carrier of weighing plate (2)
- Remove the two fastening screws (3)


- Remove fastening screw (5) from the lower part of the housing (4) and detach the upper part of the housing carefully

To install rechargeable battery:

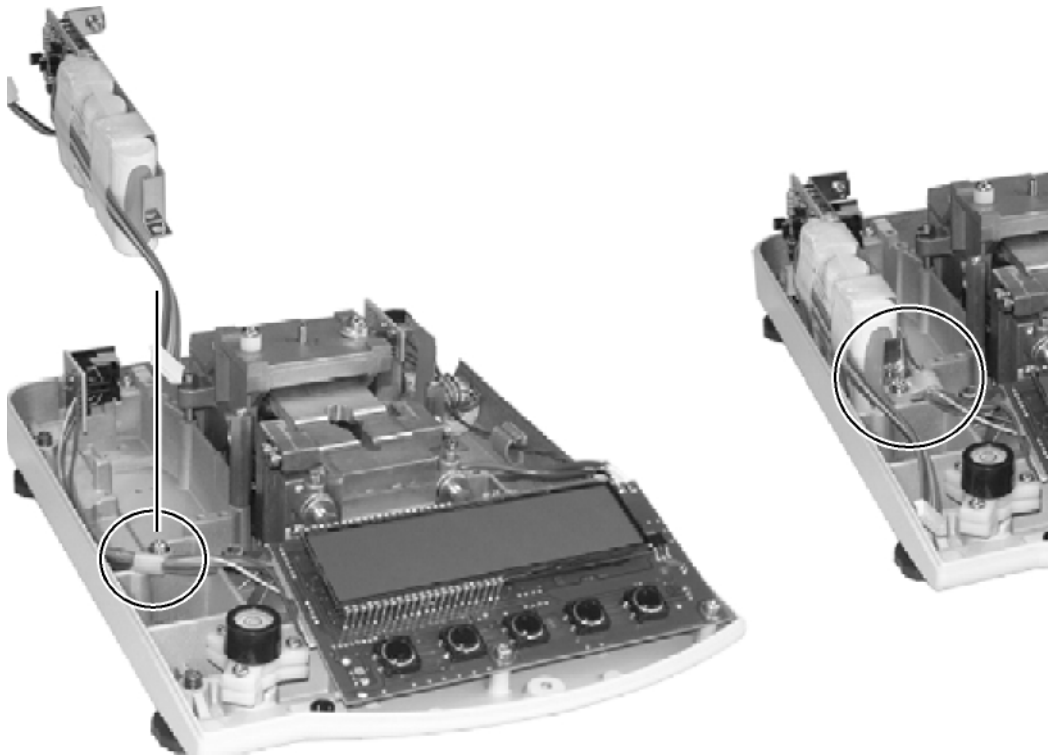- Remove the cable clip.

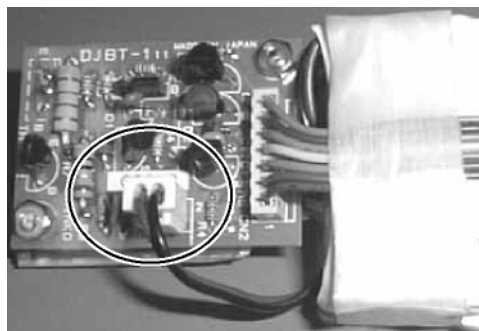
- Remove the rechargeable battery from its package and **begin by connecting the electric power supply to the circuit board of the rechargeable battery.**

- The package of the rechargeable battery is placed in the housing in a way that it can be screwed to the balance through the existing bracket by means of the cable clip loosened before.

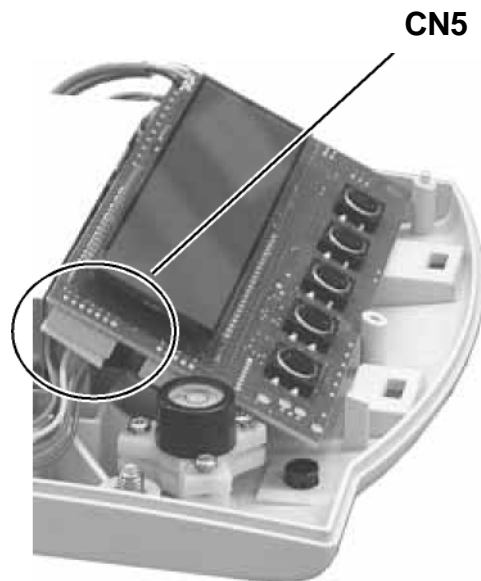- Remove the 2 fastening screws of the display circuit board
- Subsequently set up the plug-in connection to the display circuit board (3) of the balance (CN5).

- Now fasten also the display circuit board using the screws loosened previously.
- Place the upper housing section and attach it to the lower housing section by screws. Screw down both fastening screws (3) re-secure the bracket (2). Attach the weighing plate (1).

Information:

Though the rechargeable battery is immediately operative, it should be charged for at least 8 hours using the mains adaptor before being used for the first time.

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