

KERN[®] **KERN & Sohn GmbH**

Ziegelei 1
72336 Balingen-Frommern
Germany

☎ +0049-[0]7433-9933-0
FAX +0049-[0]7433-9933-149
✉ info@kern-sohn.com

Installation Instructions Interface Adapter with Cable for Blue- tooth

KERN KUP-06

Type TYKUP-06-A
Version 1.2
2023-06
GB



The current version of these instructions can also be found online under:
<https://www.kern-sohn.com/shop/de/DOWNLOADS/>
under the rubric Instruction manuals

TYKUP-06-A-IA-e-2312



Interface Adapter with Cable for Blue-tooth
Version 1.2 2023-06
Installation Instructions

Contents

1	Declaration of Conformity	2
2	Scope of delivery.....	2
3	General information	2
4	Installation	3

1 Declaration of Conformity

The valid Declaration of Conformity EC/UE is available at:

www.kern-sohn.com/ce

2 Scope of delivery

- Interface Adapter with Cable for Bluetooth
- Installation Instructions

3 General information

Transmit frequency: 2402-2480MHz

Maximum transmit power: < 20dBm

Cable length: 0.15 m



- Weighing data can be transferred via Bluetooth.
- **Only KERN KUP-adapters may be connected to the 15-pol-sub-D-connection of the balance!**

4 Installation

1. Switch off the appliance
2. Plug in the interface adapter at the KUP connection
3. Switch on the appliance. The interface adapter is automatically recognized by the appliance.

The KUP adapter is compatible with Bluetooth Low Energy (BLE) and is visible under "KERN BLE" with a unique MAC address for Bluetooth Master devices.

To access this, please use an appropriate software program / app which supports Bluetooth Low Energy (BLE). Applications exclusively using Bluetooth Classic (BLC) will not work.

The following profile must be adjusted:

Service UUID
0000fff0-0000-1000-8000-00805f9b34fb

Read characteristic UUID
0000fff1-0000-1000-8000-00805f9b34fb

Write characteristic UUID
0000fff2-0000-1000-8000-00805f9b34fb