

Ziegelei 1 72336 Balingen-Frommern Germany

+0049-[0]7433-9933-0

- +0049-[0]7433-9933-149
- info@kern-sohn.com

## Installation Instructions Interface Adapter with Cable for Ethernet

# **KERN KUP-04**

Type TYKUP-04-A Version 1.2 2023-06 GB





The current version of these instructions can also be found online under: https://www.kern-sohn.com/shop/de/DOWNLOADS/ under the rubric Instruction manuals

**TYKUP-04-A-IA-e-2312** 



#### Interface Adapter with Cable for Ethernet Version 1.2 2023-06 Installation Instructions

#### Contents

1	Scope of delivery	2
2	General information	2
3	Installation	2

#### 1 Scope of delivery

- Interface Adapter with Cable for Ethernet
- Installation Instructions

### 2 General information

Cable length: 0.15 m

1

- Via the Ethernet interface, weighing data can be transferred via a network.
- Only KERN KUP-adapters may be connected to the 15-pol-sub-D-connection of the balance!

#### 3 Installation

- 1. Switch off the appliance
- 2. Plug in the interface adapter at the KUP connection
- **3.** Switch on the appliance. The interface adapter is automatically recognized by the appliance.

In standard configuration the KUP adapter uses the DHCP log to intrude into the network. For inquiry of the configuration allocated via DHCP, as well as to the specific/static configuration of the IP-address, subnet-mask or the gateway, the KCP-commands **JNEx** can be used. A detailed description you will find in the "KERN Communications Protocol" manual, available in the download area on our KERN homepage (www.kern-sohn.com).

To activate KCP please observe the menu overview of your balance's operating instructions.