Vidrio Industrial POBEL, S.A.

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1.1. GENERAL INFORMATION

1. The package must be handled carefully. When you unpackage it, you should check that what it contains inside is what is indicated in the Packing List section. If any component is missing or otherwise damaged, contact the dealer immediately.
2. Please, read the instructions before installing the apparatus. The instruction manual will be an inseparable part of the product; For this reason, keep it in good condition for future reference.
3. The modification of the components of the product transgresses the directive of use 2009/104 / CE; and therefore: the manufacturer is not responsible for the damages that this brings as damage.
4. The product must be installed on a horizontal surface and never in places where there are vapors of contaminating or flammable substances.
5. Please, dispose of at least the first distillation before using the equipment.
6. Conservative packaging during the warranty period.
7. Warning and Observation are indicated in you can see below:

Warning: Must be strictly followed to avoid physical damage.

Note: Contains suggestions and restrictions for the product.

1.2. PACKING LIST

Description	Qty.
Metallic water still	1
Protector de salida CO ₂	1
Feed water tuve	1
Filter	1
Silicone tube 12x18	1
Silicone tube 8x15	1
Manual	1

1.3. PRODUCT SPECIFICATIONS

Specifications	10DETA0040	10DETA1040
Distillation capacity (L/H)	4	
Tank capacity (L)	10	
Cooling water (L/MIN)	1	
Conductivity 20°C (µS/cm)	1,5	
Weight (Kg)	24	
Dimensions, height x width x depth (mm)	500 x 630 x 320	
Safety	Thermostat, level switch and electrovalve	
Wattage (W)	3000	
Voltage (V)	230	110
Intensity (A)	13	27

1.4. INSTALLATION (fig. 1, 3 and 4)

1. Connect the equipment to a power outlet using the power cord CA; according to the voltage indicated in the technical specifications section.
2. Connect the grey hose to the water supply socket M. The other end of the hose, with the filter previously installed, will be connected to a $\frac{3}{4}$ "tap.
3. The drain outlet pipe D should be placed in such a position that excess water falls freely to the outside into a waste pipe about 25 cm below.
4. Place the outlet protector for CO₂ in the hole O designed for it, located on the lid of the tank A.
5. Check that tap G for cleaning is closed before and during the distillation process.

Note: the tap G is only opened for the cleaning step.

CA S LI B

fig.1

OUTLET PROTECTOR CO₂

REARME

fig.2

EB

A

O

fig.3

D

G

M

fig.4

1.5. START-UP (fig. 1 and 3)

1. Check that the installation has been carried out properly.
2. Turn the switch S to ON.
3. By opening the tap that we have connected to a ¾ "outlet, feed water. A level switch will activate when the appropriate height is reached and at that moment the green indicator L1 illuminates and remains so for the duration of the distillation.
4. Distilled water is collected in tank A and when it is filled, a level switch detects it by activating a solenoid valve that closes the water inlet avoiding overflow; at that moment the heater is disconnected, stopping the overheating of the equipment; L1 will switch off.
5. Open the tap B in order to deliver distilled water.

1.6. MAINTENANCE (fig. 1, 3 and 4)


Warning: always disconnect the power supply before cleaning

1. With switch S in the OFF position; and the CA cable disconnected from the mains, remove the upper lid of the water still.
2. Remove the lid from the boiler EB.
3. Open tap G located on the right, draining the boiler completely. When the water stops coming out, close the tap.
4. Pour in hydrochloric acid until it comes out through the drain D. It should be acting for 15 minutes.
5. Open the tap G to drain the boiler, allowing all the liquid to flow out.
6. Rinse with plenty of water and close the tap G.
7. Repeat if necessary from step 4.
8. Carry out a distillation that must be discarded.

1.7. SPARE PARTS

Description	Symbol	Cat. Number	Qty.
Contactora 230 V	C1	YE.DCONT00011	1
Contactora 110 V (10DETA1040)	C1	YE.DCONT10001	1
Electrovalve 230 V	E	YE.DELEC00002	1
Electrovalve 110 V (10DETA1040)	E	YE.DELEC00003	1
Termostato	T	YE.DTERM00001	1
Calentador metálico 230 V	R	YE.DRDES00220	1
Calentador metálico 110 V (10DETA1040)	R	YE.DRDES00110	1
Interruptor de nivel metálico, curvado	Nd	YE.DSONI00002	1
Interruptor de nivel metálico, recto	Ne	YE.DSONI00003	1
Caja reguladora de flujo	---	YP.DCAJA00002	1
Regulador de flujo 1,0 L/M	---	YP.DRECA00001	1

1.8. WATER FLOW DIAGRAM
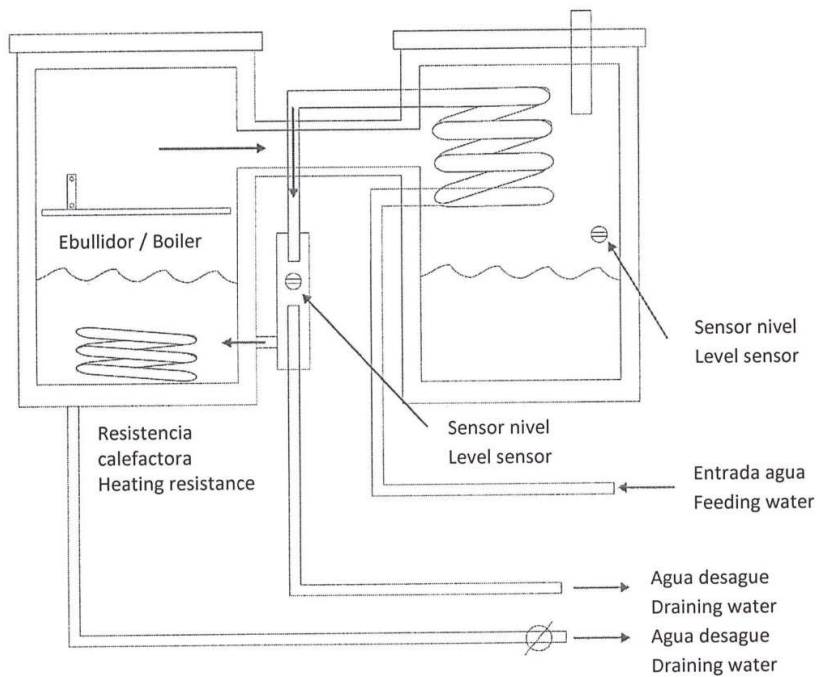


1.9. SHOOTING TROUBLE

Problem	Action
The power cord is connected to the mains, the electrical switch S is ON, and the red light on the switch does not turn on.	Measure and check that the mains voltage is 230 V (110 V for 10DETA1040) Check the status of the switch.
The water still is on, water is coming out of drain D and the green L1 light is off.	Follow this order: Remove the thermostat protector T and press inwards. Check level switches Ne and Nd. Check contactor C1. Check resistance R.
The water still is on and the water tank is not at its maximum capacity, no water circulates.	Replace the electrovalve E.
If the water level in the boiler is below the temperature sensor and the sensor triggers	Press the button REARME. (see fig.2.)

Note: It will not be allowed to repair any water still that is not properly cleaned.

1.10. ELECTRIC DIAGRAM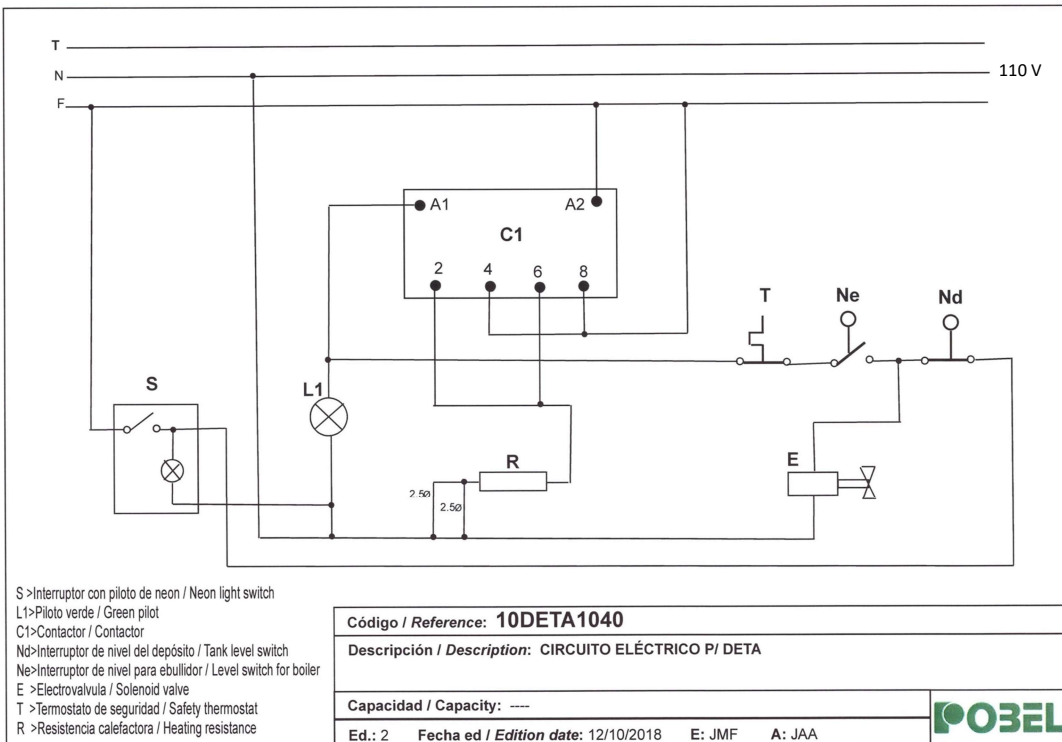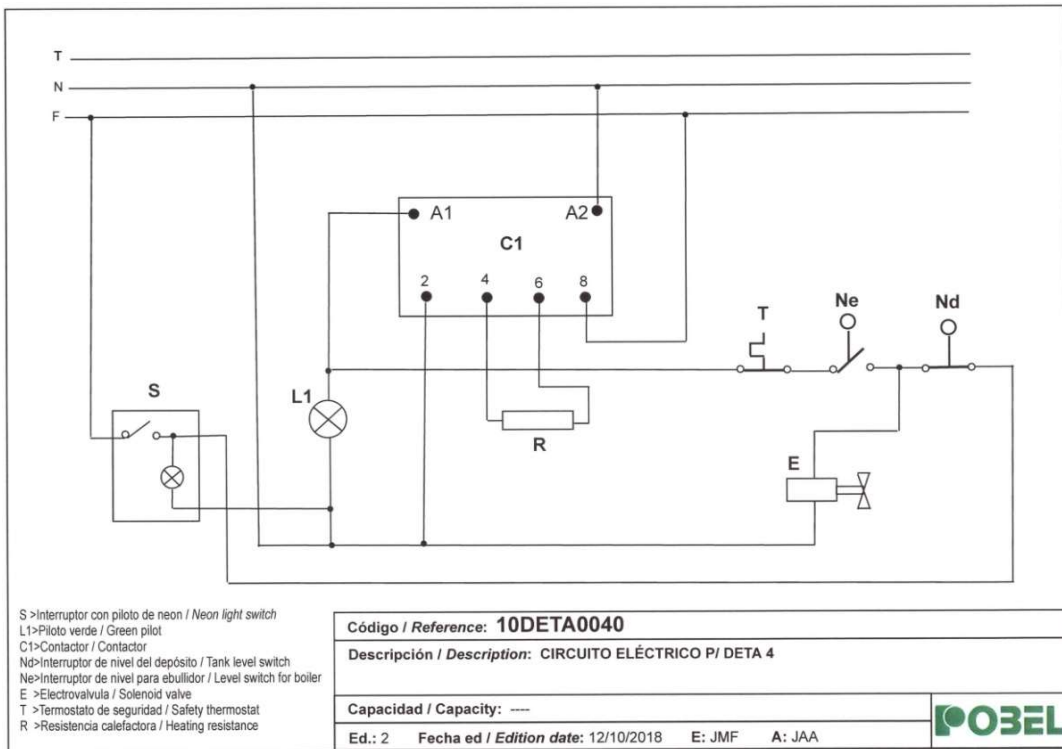

1.11. WARRANTY

1. The product covered by this Certificate is protected for two (2) YEARS after the date of the purchase delivery note of the product.
2. During this period of time, in case of manufacturing fault of the product, the user has the right to its servicing, its replacement (except if one of this options was impossible or disproportionate), to have a reduction in the price and to the resolution of the contract (when the purchaser could not require the servicing or the replacement). The servicing will be free of charge for the user. All the servicing and replacements will be carried out in a reasonable period of time.
3. The breakdowns, flaws or caused damages in the product will be excluded as consequence of: abuses, improper use, exposure to adverse weather conditions, spilled batteries, blows, falls, bad or inadequate installation, power failure, waste by normal use, servicing carried out by different people indicated by us, as well as modification of this Certificate and / or another external causes to the product.
4. The glass parts of our distillers are excluded of this warranty.
5. This warranty will be automatically cancelled in case of wrong handling of the product or there is a breach of the manual or instructions regarding the use, assemble and operation.
6. The warranty does not affect to the rights of the customer according to the law.
7. In case the user needs technical help regarding the guaranteed product, the user will contact to the distributor.
8. With the exception of choice of law rules under Spanish private international law and the provisions of the UN Convention on Contracts for the International Sale of Goods ("UN-CISG"), the substantive law of Spain shall apply to all legal relationships and transactions established by this purchase agreement.
9. Please, should give copy of the delivery note or purchase invoice with the Warranty.